



Caron Products & Services
OPERATIONS MANUAL



INSECT & ANIMAL REARING CHAMBERS

*Models: 7340-25 / 7340-33
7350-25 / 7350-33*

Dear Valued Customer:

Thank you for purchasing CARON Products & Services equipment. We appreciate your business and look forward to being your preferred supplier of controlled environment equipment products in the future.

At CARON, we are committed to continuous quality improvement. Our goal is to supply our customers with highly reliable equipment at a fair price. In order to openly monitor our performance, we would appreciate your feedback on our products and services.

If you have questions, or any suggestions for improvement based on the installation or operation of the equipment you have purchased, please contact our service department at www.caronproducts.com or 740-373-6809.

Thanks again for your business!

Revision Log

| Version | Date | Description |
|---------|----------|--|
| Rev A | 1-9-18 | Initial Release |
| Rev B | 02-12-18 | Added HUMD311-1,-4 option |
| Rev C | 04-04-18 | Edited troubleshooting section, port stoppers |
| Rev D | 07-17-18 | Updated PM Kit info |
| Rev E | 05-14-19 | Added analog output adjustable temperature range feature. |
| Rev F | 1-15-20 | Updated for flammable refrigerant |
| Rev G | 01-21-21 | Added adjustable alarm delay |
| Rev H | 03-26-21 | Changed warranty to 2 years |
| Rev J | 03-31-21 | Updated Declaration of Conformity document |
| Rev K | 04-13-21 | Updated Declaration of Conformity document |
| Rev L | 04-21-21 | Added DLOG301and RAMP301 option |
| Rev M | 11-03-21 | Updated pictures of Analog Outputs and Remote Alarm Contacts |
| Rev N | 12-30-21 | Added UKCA Declaration of Conformity |

TABLE OF CONTENTS

| | |
|--|-----------|
| Warranty | 6 |
| Equipment Overview | 10 |
| Installation..... | 13 |
| Unpacking | |
| Choosing a Location | |
| Preliminary Cleaning | |
| Installing the Port Stopper | |
| Installing the Shelves | |
| Leveling the Unit | |
| Connecting the Drain Line | |
| Connecting the Water Supply (HUMD311-1,-4 option) | |
| Connecting Electrical Power | |
| Optional Accessory Installation | 20 |
| Connecting Alarm Contacts (ALRM302) | |
| Connecting Analog Outputs (OUTP301, OUTP302) | |
| Installing Drain Water Pump (PUMP301) | |
| Operation..... | 23 |
| Using the Keypad | |
| Learning the Touch Screen | |
| Changing the Temperature Setpoint | |
| Changing the Humidity Setpoint | |
| Trend Graph | |
| Adjusting Fresh Air Ports (PORT305) | |
| Optional Accessory Operation | 31 |
| Operation of the Data Logging System. (DLOG301) | |
| Operation of the Humidity System. (HUMD311-1,-4) | |
| Operation of the LED Lighting (LGHT324, LGHT340, LGHT341, LGHT342) | |
| Operation of the Front 6" Mounted Recorders (RCDR316, RCDR317) | |
| Operation of the Front 10" Mounted Recorders (RCDR318, RCDR319) | |
| H2O2 sterilization cycle (STER305) | |
| Operation of the Ramp and Soak Cycle (RAMP301) | |
| Calibration | 62 |
| Calibrating the Temperature | |
| Calibrating the Humidity | |
| Calibrating Optional Chart Recorders | |
| Alarms | 66 |
| Alarm System Overview | |

| | |
|---|------------|
| Snoozing the Speaker | |
| Muting the Speaker | |
| Changing Alarm Setpoints and Delay | |
| Alerts | 72 |
| Alert System Overview | |
| Maintenance | |
| Info | 75 |
| Info System Overview | |
| Advanced Features | 76 |
| Setting the time & day | |
| Door Heat | |
| Humidity Control | |
| Locking the controls | |
| Changing the passcode | |
| Factory menu & troubleshooting | |
| Preventative Maintenance | 92 |
| Specifications | 94 |
| Electrical Schematics | 95 |
| Troubleshooting | 104 |
| Spare Replacement Parts | 105 |
| Appendix A – Declaration of Conformity | 107 |

WARRANTY INFORMATION

EQUIPMENT LIMITED WARRANTY

Please review this section before requesting warranty service. At CARON, one of our primary goals is to provide customers with high levels of personal service and top quality products, delivered on time, backed by technical service and supported for the life of the product.

Before contacting us for warranty service, please be aware that there are repairs that are not covered under warranty.

WARRANTY DEFINED

Caron Products & Services, Inc. (herein after CARON) hereby warrants that equipment manufactured by CARON is free from defects in materials and workmanship when the equipment is used under normal operating conditions in accordance with the instructions provided by CARON.

COVERED:

- Parts and labor for a period of two (2) year from date of shipment.
- Any part found defective will be either repaired or replaced at CARON's discretion, free of charge, by CARON in Marietta, OH. Parts that are replaced will become the property of CARON.
- If CARON factory service personnel determine that the customer's unit requires further service CARON may, at its sole discretion, provide a service technician to correct the problem, or require the return of the equipment to the factory or authorized service depot.
- CARON will have the right to inspect the equipment and determine the repairs or replacement parts necessary. The customer will be notified, within a reasonable time after inspection, of any costs incurred that are not covered by this warranty prior to initiation of any such repairs.

NOT COVERED:

- Calibration of control parameters.
- Improper installation; including electrical service, gas and water supply tubing, gas supplies, room ventilation, unit leveling, facility structural inadequacies or ambient conditions that are out of specification.
- Cost of express shipment of equipment or parts.
- Any customer modifications of this equipment, or any repairs undertaken without the prior written consent of CARON, will render this limited warranty void.
- CARON is not responsible for consequential, incidental or special damages; whether shipping damage or damages that may occur during transfer to the customer's point of use. When the equipment is signed for at the customer's site, ownership is transferred to the customer. Any damage claims against the shipping company become the responsibility of the customer.
- Repairs necessary because of the equipment being used under other than normal operating conditions or for other than its intended use.
- Repair due to the customer's failure to follow normal maintenance instructions.
- Parts considered consumable; including: light bulbs, filters, gases, etc.
- Damage from use of improper water quality.
- Damage from chemicals or cleaning agents detrimental to equipment materials.
- Force Majeure or Acts of God.

This writing is a final and complete integration of the agreement between CARON and the customer. CARON makes no other warranties, express or implied, of merchantability, fitness for a particular purpose or otherwise, with respect to the goods sold under this agreement. This warranty cannot be altered unless CARON agrees to an alteration in writing and expressly stated herein shall be recognized to vary or modify this contract.

Ohio Law governs this warranty.

EQUIPMENT INTERNATIONAL LIMITED WARRANTY

Please review this section before requesting warranty service. At CARON, one of our primary goals is to provide customers with high levels of personal service and top quality products, delivered on time, backed by technical service and supported for the life of the product.

Before contacting your distributor for warranty service, please be aware that there are repairs that are not covered under warranty.

WARRANTY DEFINED

Caron Products & Services, Inc. (herein after CARON) hereby warrants that equipment manufactured by CARON is free from defects in materials and workmanship when the equipment is used under normal operating conditions in accordance with the instructions provided by CARON.

COVERED:

- Parts for a period of two (2) years from date of shipment.
- Any part found defective will be either repaired or replaced at CARON's or their authorized representative's discretion. Parts that are replaced will become the property of CARON.
- If CARON or their authorized representatives determine that the customer's unit requires further service, CARON or the representative may, at its sole discretion, provide a service technician to correct the problem, or require the return of the equipment to the an authorized service depot.
- CARON or their authorized representative will have the right to inspect the equipment and determine the repairs or replacement parts necessary. The customer will be notified, within a reasonable time after inspection, of any costs incurred that are not covered by this warranty prior to initiation of any such repairs.

NOT COVERED:

- Calibration of control parameters.
- Improper installation; including electrical service, gas and water supply tubing, gas supplies, room ventilation, unit leveling, facility structural inadequacies or ambient conditions that are out of specification.
- Cost of express shipment of equipment or parts.
- Any customer modifications of this equipment, or any repairs undertaken without the prior written consent of CARON, will render this limited warranty void.
- CARON and their representative are not responsible for consequential, incidental or special damages; whether shipping damage or damages that may occur during transfer to the customer's point of use. When the equipment is signed for at the customer's site, ownership is transferred to the customer. Any damage claims against the shipping company become the responsibility of the customer.
- Repairs necessary because of the equipment being used under other than normal operating conditions or for other than its intended use.
- Repair due to the customer's failure to follow normal maintenance instructions.
- Parts considered consumable; including: light bulbs, filters, gases, etc.
- Damage from use of improper water quality.
- Damage from chemicals or cleaning agents detrimental to equipment materials.
- Force Majeure or Acts of God.

This writing is a final and complete integration of the agreement between CARON and the customer. CARON makes no other warranties, express or implied, of merchantability, fitness for a particular purpose or otherwise, with respect to the goods sold under this agreement. This warranty cannot be altered unless CARON agrees to an alteration in writing and expressly stated herein shall be recognized to vary or modify this contract.

Ohio Law governs this warranty.

Caron Products & Services, Inc.
PO Box 715 · Marietta, OH 45750
740-373-6809

INTERNATIONAL SYMBOLS AND DEFINITIONS



Help



Information



Warning of hazardous area



Warning of hot surface



Warning of dangerous electric voltage



Warning of risk of fire



Earth (ground) protective
conductor

WARNINGS



Local government may require proper disposal

TO LOCATE REFRIGERANT TYPE AND PRESSURES, SEE SERIAL LABEL LOCATED ON THE OUTSIDE OF THE UNIT

FOR HYDROCARBON (R290 PROPANE) REFRIGERANT UNITS

R290 is highly flammable and must be treated with proper care.



Do not damage the refrigeration circuit. Do not store explosive substances in the unit. Component parts shall be replaced with like components and servicing shall be done by authorized personnel to reduce the risk of possible ignition.



DANGER – Flammable Refrigerant Used. Risk of fire or explosion.

- Do not puncture refrigerant tubing
- Do not use mechanical devices to defrost refrigeration equipment
- Unit to be repaired only by trained service personnel



CAUTION – Flammable Refrigerant Used. Risk of fire or explosion.

- Consult repair manual, owners guide before attempting to service this product. All safety instructions must be followed.
- Dispose of properly in accordance with federal or local regulations.



CAUTION – Do not use any electrical appliance within the Environmental Chamber, other than those recommended by the manufacturer.

EQUIPMENT OVERVIEW

Congratulations! You have just purchased the latest technology in Insect/Animal Rearing chambers. Before using the equipment, familiarize yourself with key components of the product and thoroughly read this manual.

This user's manual covers both 7340 Insect Rearing and 7350 Animal Rearing Series Chambers




Model 7340




Model 7350

EQUIPMENT OVERVIEW – CONTINUED

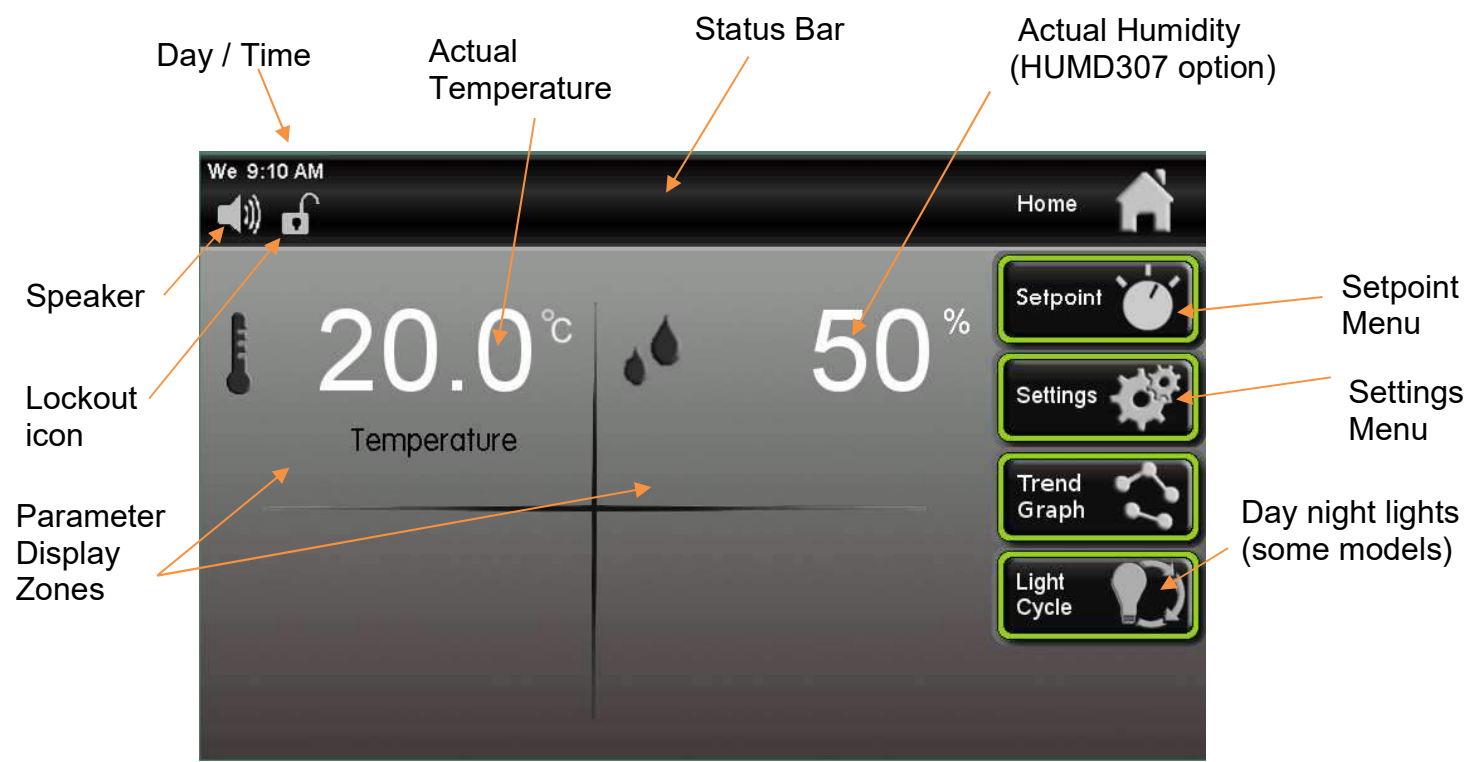

(Home Screen)

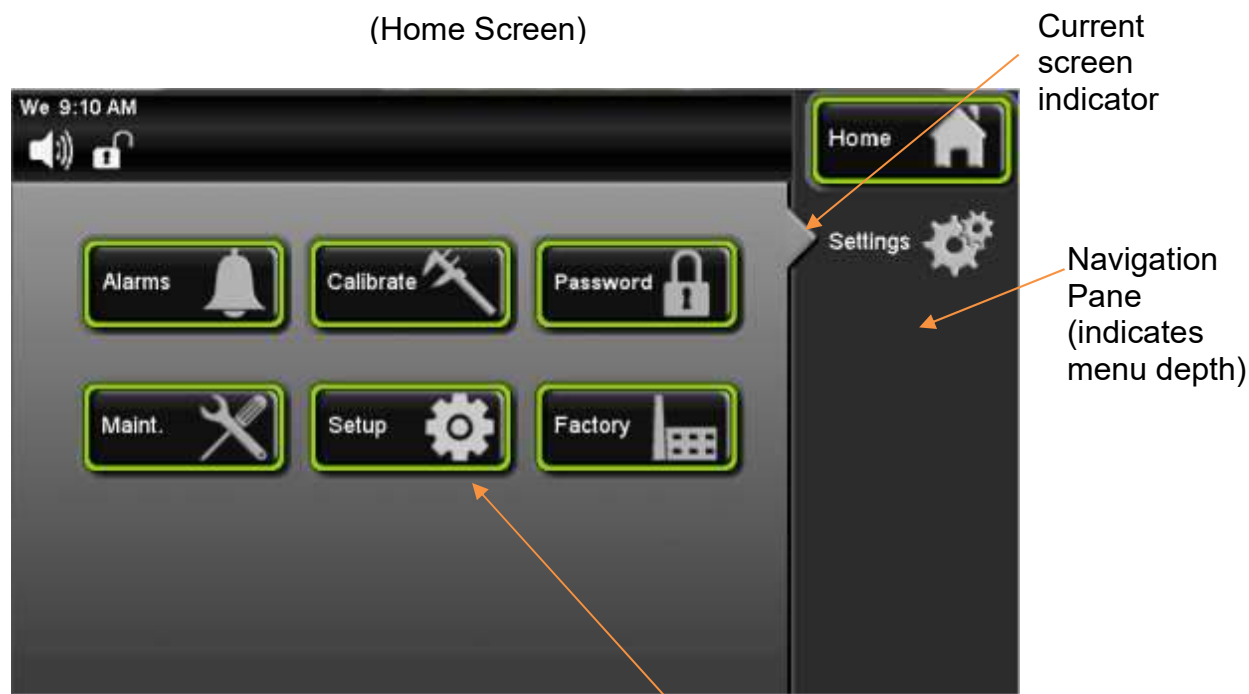

(Settings Screen)

INSTALLATION

Unpacking

Your new unit has been thoroughly packaged to avoid shipping damage. However, the unit should be fully inspected upon arrival before signing for receipt. If the package has visual damage, notes should be made on the freight bill and signed by the delivery company. In the event of concealed damage after the unit is uncrated, keep the carton and packaging material. Call the shipping company within 7 days of receipt, request inspection and retain a copy of the inspection report.

For prolonged periods of inactivity leave the unit unplugged and securely stored.

Caron provides full on-site installation services for all models. Our installation services guarantees the proper set-up and startup of all equipment. Please contact the Service Department at 740-373-6809 or www.caronproducts.com for details.

For detailed instructions on how to safely remove the chamber off the shipping pallet, see document located on the chamber.

Choosing a Location



This product weighs in excess of 500 pounds. Ensure that sufficient resources are available to safely move the product.

To ensure proper operation, the unit must be located on a firm level surface, capable of supporting approximately 900 pounds. The unit should be located in an 18°C to 25°C ambient area and where there is no direct airflow from heating and cooling ducts as well as out of direct sunlight. Allow four inches of clearance on all sides of the product to allow for connections and airflow. The unit is designed to be used under the following conditions:

- Indoor use only
- Altitude up to 2000m.
- Maximum relative humidity: non-condensing
- Mains supply voltage fluctuations up to +/- 10% of the nominal voltage; damage may occur if voltage varies more than 10%
- Transient overvoltages up to the levels of overvoltage category II
- Temporary overvoltages occurring on the mains power supply
- Pollution degree: 2
- Ingress protection: IPXO

* Depending on user setpoints, these units may not need a drain. Drains are recommended at temperatures below 15°C.

The unit requires a dedicated electrical connection. Power requirements vary depending upon the chamber model, see Connecting Electrical Power section.

Choose a location where these facilities are, or can be made available. If a water source or a drain is not available, contact CARON customer service and ask about our CRY5102 product line or click this web link for information on the product:

www.caronproducts.com

Preliminary Cleaning

Your new environmental chamber was thoroughly cleaned prior to leaving the factory. It is recommended, however, to clean all interior surfaces with a general purpose laboratory cleaning agent to remove any shipping dust or dirt prior to using the product. Contact Caron if there is any doubt to the compatibility of the cleaning agent being used with the chamber. After cleaning, dry all interior components with a sterile cloth as necessary.

Installing the Port Stoppers

The unit has an access port built into each side of the cabinet. The ports are designed to allow customer access for equipment validation and for installation of other equipment inside the chamber. These ports should be sealed with the provided rubber stoppers to allow the chamber to function properly. Install the stoppers provided in the port on each side of the unit.



R290 REFRIGERANT UNITS

DANGER – Flammable Refrigerant Used. Risk of fire or explosion.

- No equipment that uses an open flame should be placed inside the unit.
- Do not use instrumentation or equipment that incorporates potential ignition sources, e.g. open contact switching, brushed DC and AC motors, etc.
- Do not use electrical appliances within the unit, other than those recommended by the manufacturer.

Installing the Shelves

Each new 7340-25/7350-25 Series chamber includes 4 perforated stainless steel shelves. Each new 7340-33/ 7350-33 Series environmental chamber includes five perforated stainless steel shelves. Each shelf requires two shelf channels for installation. The left and right shelf tracks are the same. Prior to installation, take time to consider what the size of the product being placed in the chamber will be and set the shelf spacing accordingly. Additional shelving can be purchased through CARON customer service if necessary. The chamber should be empty when being moved.

To install the shelf channels insert the rear tab on the shelf channel into the rear wall on the side wall of the chamber. Then insert the front tab into the front pilaster. Push the entire shelf channel towards the rear of the unit and snap it down into place.



Each shelf is capable of supporting a uniformly distributed load of 50 pounds. The maximum chamber capacity (stationary) is: 500 pounds.



Do not have multiple loaded shelves out simultaneously or the chamber may tip.

Leveling the Unit

Place a level on the middle shelf of the incubator. Adjust the leveling feet until the unit sits level left to right and front to back. Even if the unit is level without adjustment, the leveling feet should still be lowered to avoid the cabinet moving while opening and closing the outer door & prevent a flat spot from forming on the casters.



Connecting the Drain Line



When using a pressurized water source, failure to connect the unit to a drain could result in facility flooding.



The chamber drain connection is located in the bottom middle of the back of the chamber. A 3/8" tube fitting, tubing and wire ties are supplied in the unit parts kit. Insert the tube fitting into the tubing, secure tubing to fitting with provided wire tie. Insert fitting into drain connection. Pull on the tubing after installation to make sure it is secure. Route the drain tubing to a local floor drain. Duplicate fitting installation on other end of tubing if necessary



The drain line relies on gravity to remove water from the chamber. The drain line must remain below the chamber to drain properly. Kinks or elevations in the drain line above the cabinet drain will not allow the chamber to drain.

If a local floor drain is not available, a variety of accessories are available through CARON customer service. These accessories can also be viewed at www.caronproducts.com.

For non-humidified chambers operated above 15°C, a drain is not needed. Install the 3/8" plug into the drain connection.

Connecting the Water Supply (HUMD311-1, -4 option)

This applies to units that have the HUMD311-1,-4 humidity option.


To ensure proper operation, distilled or deionized water is required as a supply on units that have humidity control. If these water sources are not available contact CARON customer service.



Use only distilled or deionized water with a resistivity between 50K Ω -CM and 1M Ω -CM and a pH of greater than 6.5. Using water outside this range will void your warranty.



Do not use water that contains chloramines. Chloramines can damage internal rubber gaskets resulting in leaks.

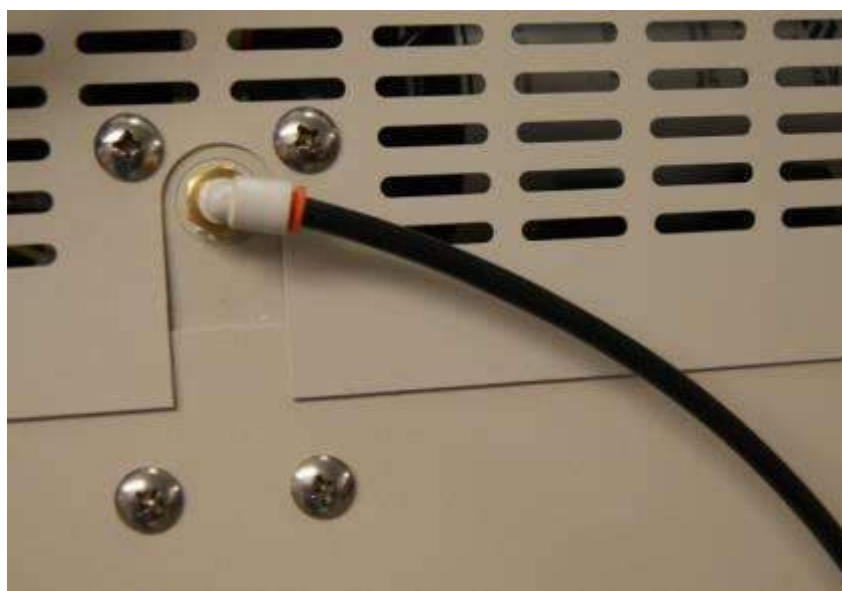

A water inlet fitting on the back of the unit and ¼" black tubing are provided to connect the water supply to the chamber. Connect an appropriate water supply to the fitting. Incoming line pressure should be regulated to not exceed 80 psi.

If a Condensate Recirculator water recycling system was purchased as a water supply, refer to its user's manual for proper installation.

Connecting Electrical Power



Connect each chamber to a grounded circuit. Failure to do so could result in electrical shock.

The unit requires a dedicated electrical outlet. See table below for model specific power required and connection.

| Model # | Power Requirements | Plug Connection |
|---------|---------------------|-----------------|
| -1 | 115V, 60Hz, 16A FLA | NEMA 5-20 |
| -2 | 230V, 60Hz, 10A FLA | NEMA 6-15 |
| -3 | 230V, 50Hz, 8A FLA | CEE 7/7 |

When the required electrical connection is available, plug the provided power cord into the unit and the electrical outlet.



The mains power supply cord must meet the requirements listed above. The use of an inadequate mains power supply cord could result in equipment failure or personal harm to the user.



In the event of a power outage the unit will automatically restart when the power is restored.


OPTIONAL ACCESSORY INSTALLATION

Connecting Alarm Contacts (ALRM302)

With the purchase of ALRM302, a set of terminals on the rear of the unit is provided to monitor temperature and humidity alarms.

With the alarm contacts, the terminals provided allow for a NO (normally open) output, a NC (normally closed) and COM (common) connection. In the event of an alarm condition or power failure, the NO contact will close, and the NC contact will open. Once the alarm is cleared, the contacts return to their normal conditions. Insert the appropriate wire into the terminal and tighten down the screw terminal on top of the connector.

| Terminal Connection | Unit off | Normal | Alarm |
|---------------------|----------|--------|--------|
| N/O to C | Closed | Open | Closed |
| N/C to C | Open | Close | Open |

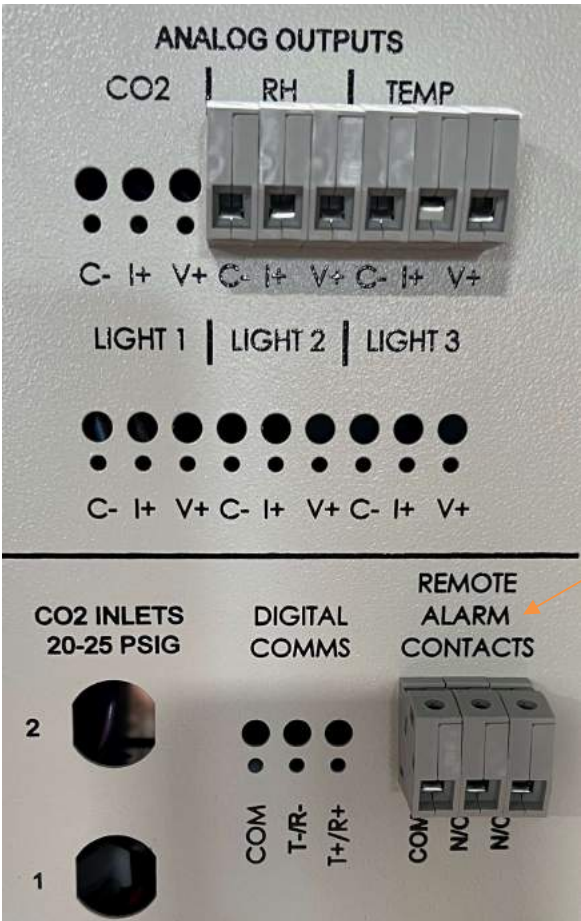


Alarm Contact Connections


Connecting Analog Outputs (OUTP301, OUTP302)

With the purchase of OUTP301, OUTP302 the controls are equipped with analog outputs. A set of terminals are provided on the back of the unit to connect to analog outputs.

Analog Output Connections

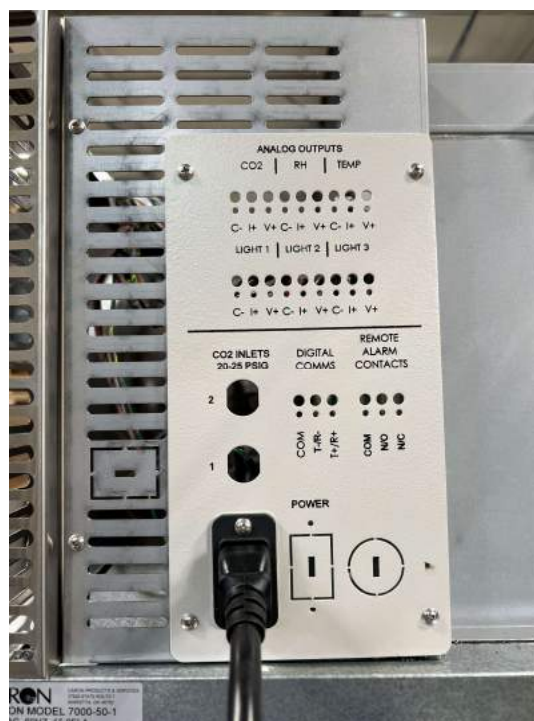

Analog outputs provide either a milliamp (4-20mA) or voltage (0-5V) signal output to represent each of the displayed temperature (and humidity) values. These options can be used for connection to in-house data acquisition, recorder, or alarm system. The temperature parameter (only) is adjustable in its scaling and is accessible at the Analog Output screen.

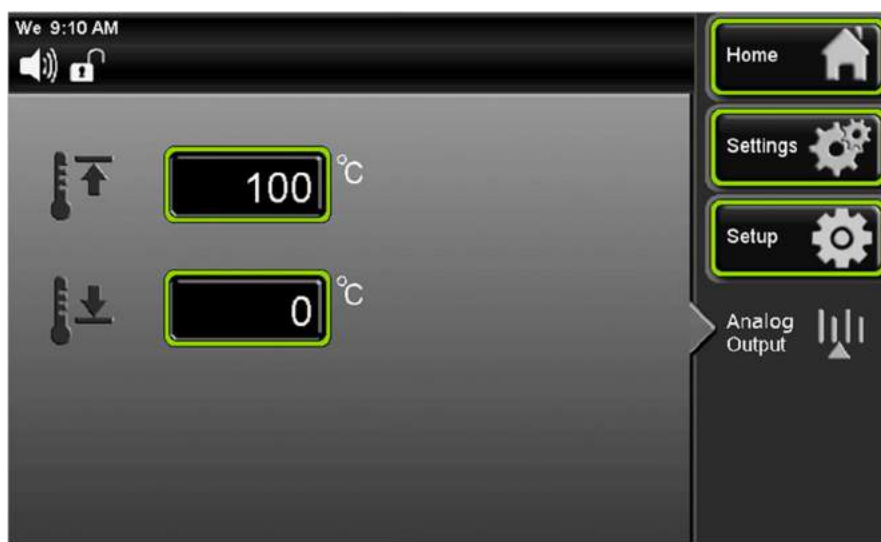


| Parameter | Voltage | Current | Corresponding Value |
|-----------------|---------|---------|----------------------------|
| Temperature | 0 – 5 V | 4-20 mA | -50 – 100 °C (adjustable*) |
| Humidity | 0 – 5 V | 4-20 mA | 0 – 100 %rh |
| CO ₂ | 0 – 5 V | 4-20 mA | 0 – 20 %CO ₂ |

*Default range is -50C to +100C. Temperature scale low range is adjustable from -50C to 0C. Temperature scale high range is adjustable from 1C to 100C.

Connect shielded wires to the appropriate signal terminals: I(+) for current (mA) or V(+) for voltage (DC). For both current and voltage outputs, COM(-) is common terminal.

Installing Drain Water Pump (PUMP301)

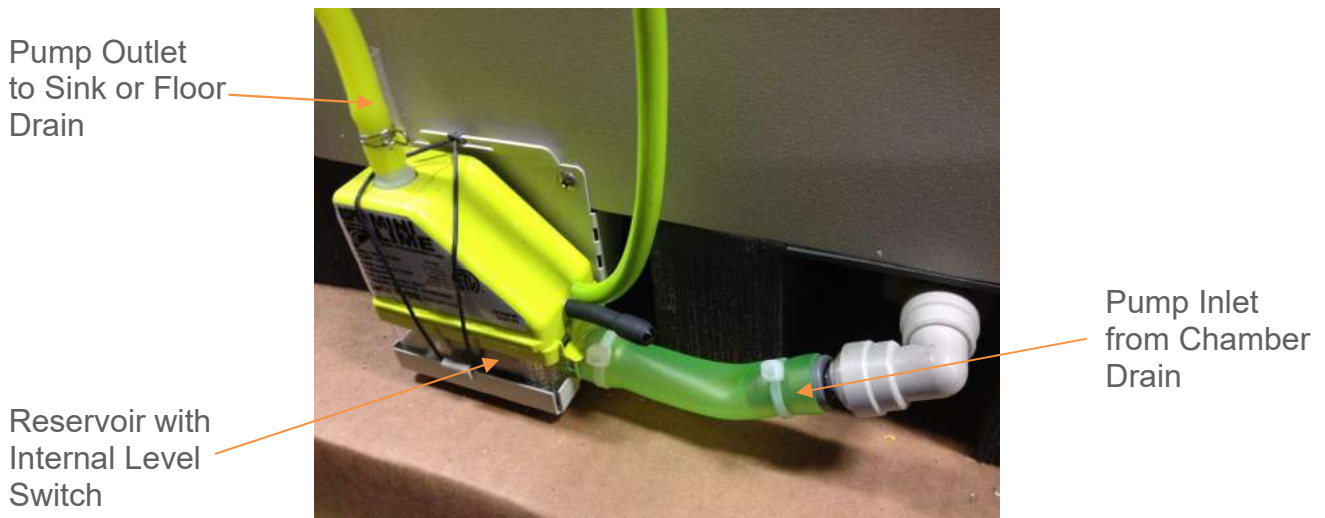

In applications where a floor drain is not available and a CARON water recycling system is not being used, a drain pump can be purchased to pump any excess condensate from the chamber to a local sink or drain. The pump is located near the middle of the back of the chamber. Connect the supplied tubing from the pump to the sink / drain. The tubing may be run vertically into a ceiling but should not exceed 15 feet height. The pump is equipped with a small reservoir on the bottom of the pump with an internal level switch that will automatically turn the pump *ON* when it is full to drain the water out of the reservoir and into a floor or sink drain.


OPERATION

Before the incubator can be commissioned for use, make sure that the following steps have been completed:

- Chamber is properly installed and level.
- The appropriate utilities connected to the chamber.


With the above mentioned steps complete, the power switch located on the right side, near the top of the unit exterior, can be turned on.

Within a few minutes, the temperature will begin to approach setpoints. Here is an overview of the home screen.

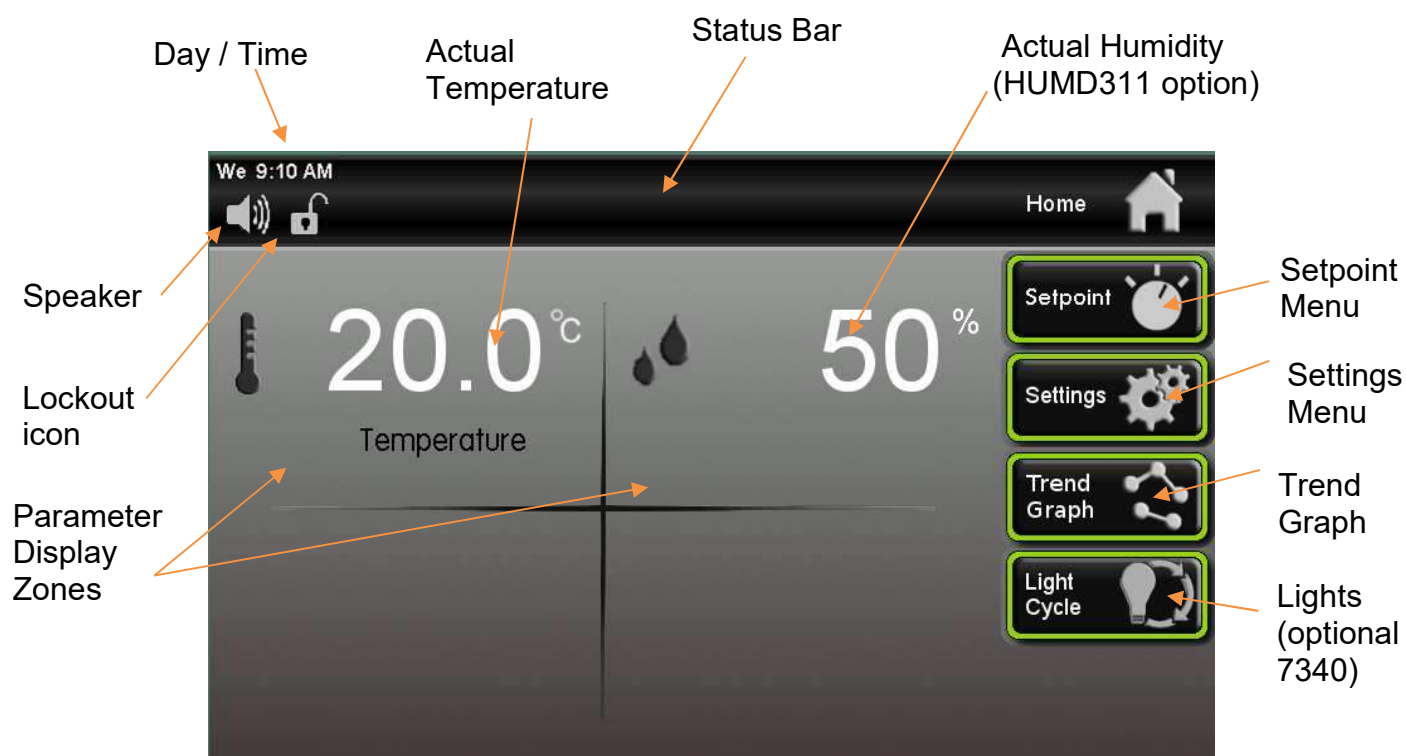

Using the Keypad

This control system uses a numeric keypad to enter all parameter values. Similar to a calculator, this allows quick and precise entry of values. When any numeric value button is pressed, the keypad display will pop up over the current display.

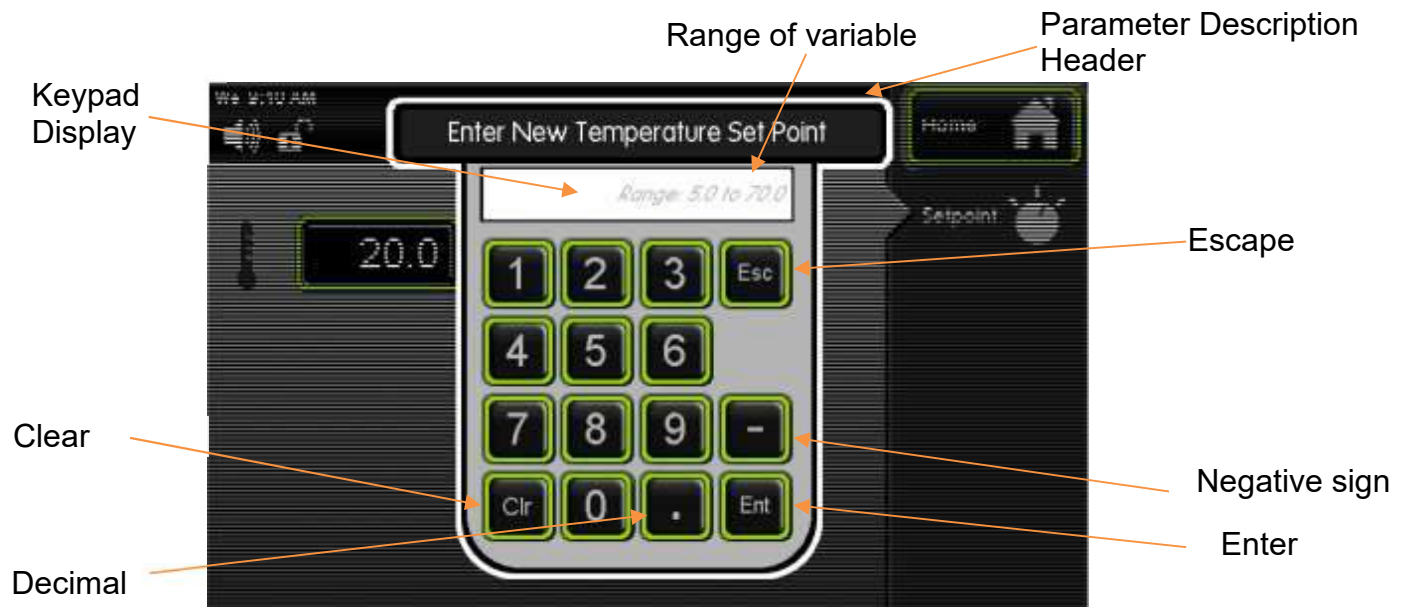

The Parameter Description Header tells what parameter is being changed. The Keypad Display shows allowable values of the parameter being changed (initially) and displays the entered value (when a button is pressed).

The Escape “Esc” button aborts the entry and returns to the previous screen without changing the value. The Clear “Clr” button erases the value that you have entered. After you have entered the value that you want, pressing the Enter “Ent” button and the new value will take affect. This also closes the keypad window. Other keypad buttons include a decimal point button and negative button.

If an invalid numeric button is pressed such that it would create an entry above the parameter’s range, the entered number will not display. For example, if the temperature setpoint range is 5.0 to 70.0, pressing ‘8’ followed by an ‘0’, only the ‘8’ will display.


If an invalid entry is made with an entry below the range (such as a ‘4’ followed by the ‘Ent’ button), then the entry will clear and the range will be re-displayed.

Learning the Touchscreen


To save power and ensure long product life, the touchscreen display has a few features that can be changed to reduce screen brightness and initiate a Screen Saver mode.



High / Low button: high or low screen brightness, preset values.

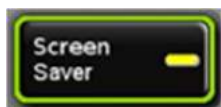



Screen Saver: By pressing the Screen Saver button “on” this will automatically enter screen saver mode after 15 minutes. At this time, the screen will be completely blank (ie. black). The illuminated Caron logo (see Equipment Overview section) shows that the unit is powered on and functioning. To wake-up the touchscreen, simply press anywhere on the touchscreen and the main screen will display. If the unit has an alarm condition, the touchscreen will not go into screen saver mode. If an alarm condition occurs while in screen saver mode, the display will automatically wake up and display the alarm.

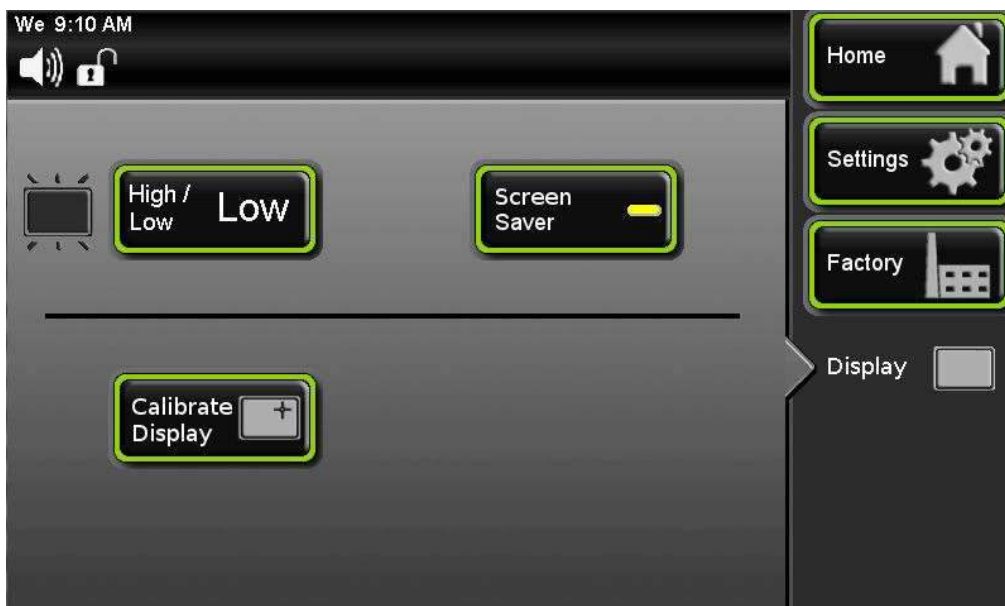



Changing the Temperature Setpoint

The steps below walk through an example of changing the temperature setpoint from 30.0 °C to 20.0 °C. This example shows humidity control as well (HUMD307 option). Here is the display of the home screen.

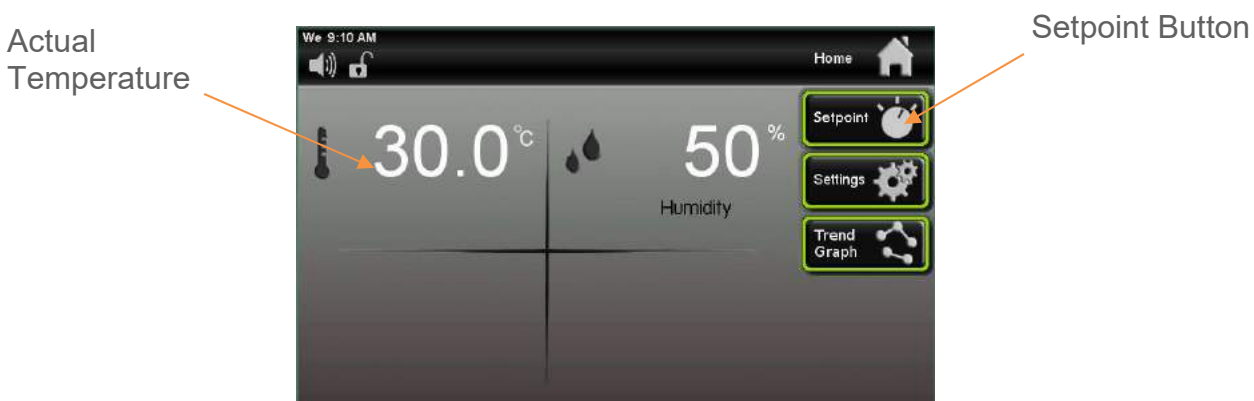

To set the temperature setpoint, press the  (Setpoint) button on the right side of the screen.

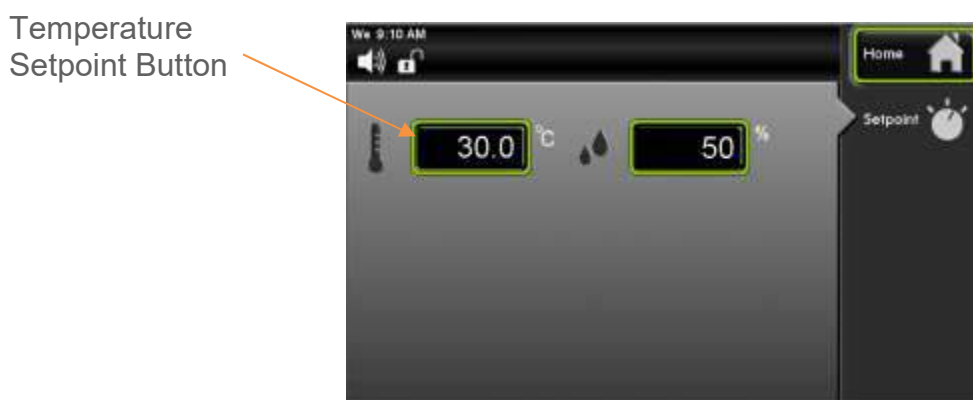


Once the Setpoint screen appears, press the  (Temperature Setpoint) button. (In this example the temperature setpoint initially has a value of '30.0'; this will vary with different initial setpoint values.)

A temperature setpoint window will appear. Enter the temperature setpoint by using the keypad. For a setpoint of 20, press '2'), then ('0'), followed by the (Ent) button. Correct any mistakes with the (Clr) and reenter the value.

Once the Enter key has been pressed, the pop-up keypad disappears and the screen returns to the Setpoint display with the new value of 20.0 °C in the temperature set point button.

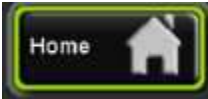

Press the  (Home) button to return to the main screen.

Changing the Humidity Setpoint (HUMD311-1,-4)


The steps below walk through an example of changing the humidity setpoint. Here is the display of the home screen.



To set the humidity setpoint, press the  (Setpoint) button on the right side of the screen




Once the setpoint screen appears, press the  (Humidity Setpoint) button.

Enter the new humidity setpoint on the keypad as desired and press  (Enter) when complete.




Home Button

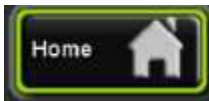

Press the (Home) button to return to the main screen.

Trend Graph

This feature allows the user to view a graph that displays the Temperature and Humidity control line over a period of time. This enables user to view the chamber's most recent performance. Trend Graph data cannot be downloaded.

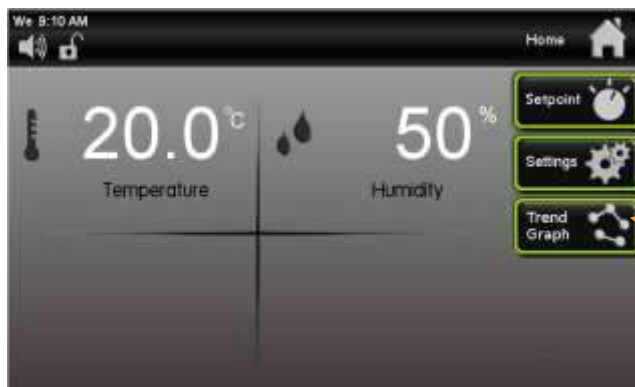

Trend Graph Button

Trend Graph displays Temperature and Humidity. Orange bar indicates current time



Adjusting the Fresh Air Ports (PORT305)

Model 7350 comes standard with these fresh air ports. The fresh air ports will allow up to 15 air exchanges per hour when fully opened. The fresh air port inlet is located at the top of the chamber, and the exhaust air port is located near the bottom of the chamber. The fresh air is drawn into the chamber plenum and distributed throughout the chamber a percentage of the air that is in the chamber exits through the exhaust port.

Both inlet and exhaust ports can be adjusted to allow variable airflow through the chamber. Both ports have a fine mesh screen to prevent objects from getting in or getting out of the chamber.



If the option STER305 sterilization cycle was purchased with model 7350 or model 7340 with the PORT305 option, make sure that these ports are closed before initiating the sterilization cycle.

ACCESSORY OPERATION



Operation of the Data Logger (DLOG301)

The DLOG301 option provides the customer with a means of logging data electronically for viewing at a later date. Logged variables are Temperature, Humidity, CO₂ and Light Intensity (but only if the chamber is equipped with those features.) All data is time-stamped with year, month, day of the month, hour, minute, 24 hour time (ISO 8601 format). This data is stored internally in the chamber in non-volatile memory.

Note: The date and time are logged within the actual file name. The file's "Date modified" field is not maintained and therefore may not reflect the actual date and time the file was created.

Data is logged every 5 minutes (provided the chamber is on); more than 10 years of data can be stored in memory. If the internal memory fills up, new data overwrites the oldest data.



Continuous writing to the flash drive necessitates a high quality industrial grade device. Use only the flash drive provided by Caron (or equivalent: single level cell memory, wear leveling algorithms, error correcting code).



File name format is "DATE START YYYY-MM-DDTHH-MM_.csv" (hours in 24 hour time)

When the chamber is on, the chamber's history data is being stored even when a flash drive is **not** inserted in the USB port. This data may be retrieved anytime using the provided USB flash drive.

Here are the methods for retrieving data:

Continuous logging of data

Insert the flash drive into the chamber's USB port. When first inserted, it creates a .csv file called 'DATA START' with the current date and time in the file name. At 5 min intervals, the chamber's process values are appended to the file. (The file will get as large as the flash drive, permitting several years of uninterrupted data storage.)




USB icon appears in in Status bar indicating that data is being written to flash drive. To retrieve the data press the 'Eject' button, then insert the flash drive into a computer to upload the data.

Upon re-insertion of the flash drive, a new .csv file is created, even if the old file is still present. File name nomenclature is "DATE START YYYY-MM-DDTHH-MM.csv".

History Retrieval





Select the 'Auto Export' feature on the USB menu screen. Insert the flash drive into the chamber's USB port. A new .csv file is automatically created on the flash drive with all the stored history data. The file name nomenclature is "DATE END YYYY-MM-DDTHH-MM.csv".

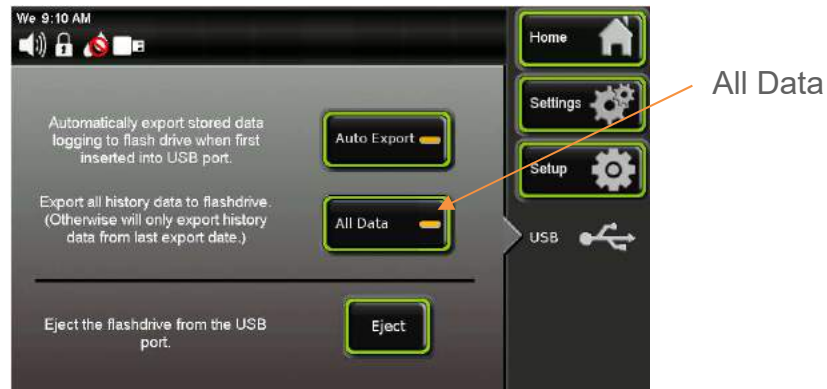

There is also an 'All Data' feature to indicate if the upload should include all data (since the unit has been used) or just the history data since a flash drive was last inserted. An 'Info' button will appear in the status bar warning the user not to remove the flash drive while the data is being uploaded. The length of time to upload the file will depend on the file size. When the 'Info' button disappears from the status bar, press the 'Eject' button to safely remove the flash drive. Now the data can be uploaded to a computer for viewing.

When using the Continuous Logging of Data method, nothing on the touchscreen has to be setup. However using the History Retrieval method will require going into the USB screen to select either the 'Auto Export' or 'All Data' buttons before inserting flash drive into USB port.

To select the 'Auto Export' and 'All Data' buttons.

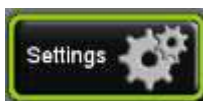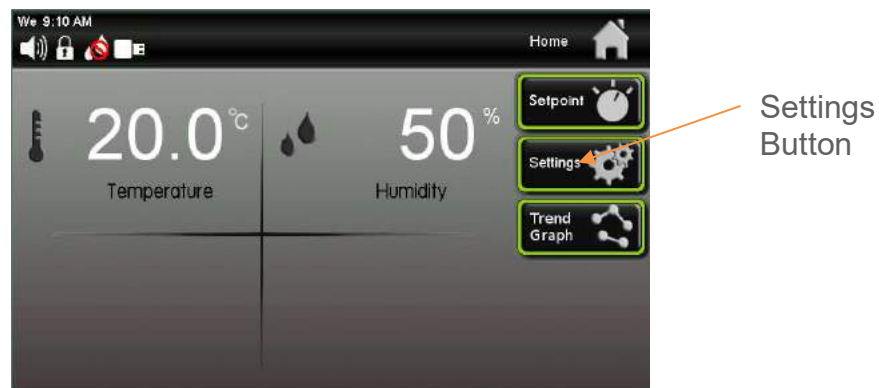

Press the (Settings) button.



Setup Button




Press the (Setup) button.



USB Button



Press the (USB) button.

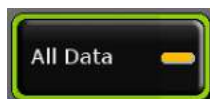


When the 'All Data' button is selected this will retrieve data starting at the point of the last download, and continuing to the present time.



When the 'Auto Export' button is selected this will retrieve the data starting at the point of the last download, and continuing to the present time.

USB flash drive icon

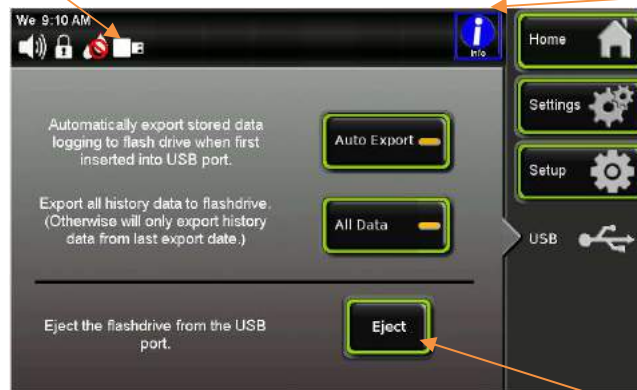

Info button

Eject button

When flash drive is inserted into the USB port a 'USB flash drive' icon and flashing 'Info' button appears in the status bar indicating that the data is being downloaded to


the flash drive. Once 'Info' icon stops flashing select the 'Eject' button.



Wait until the USB icon disappears to safely remove the flash drive from the USB port.

Note: Press the Eject button before removing the flash drive from the chamber, otherwise there could be the risk of corrupt data.

Here is a graphic to illustrate how the data retrieval works.

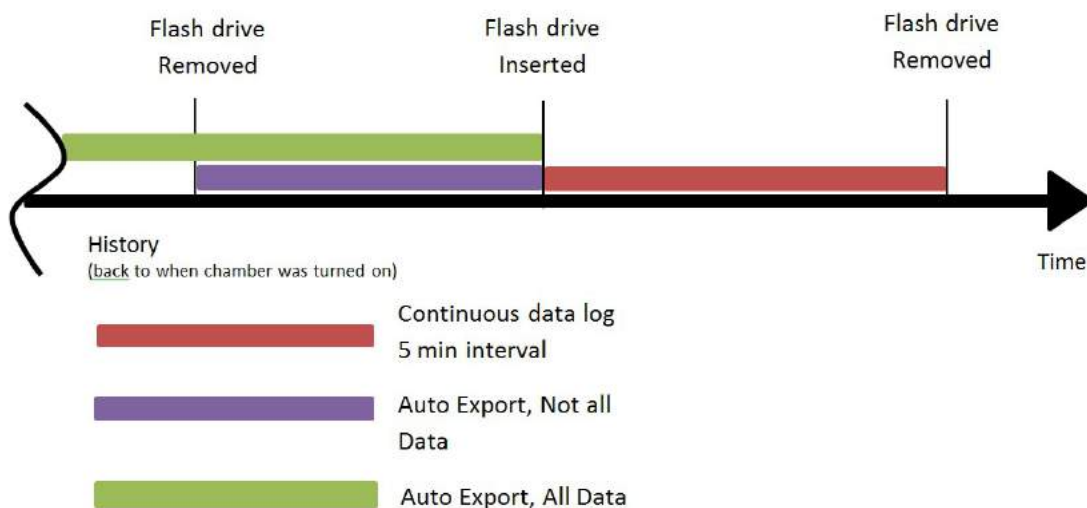

HUMD311-4,-4 Humidity System

The HUMD311 Humidity option has controlled humidity injection. Typically dehumidification is done by utilizing a wet coil (dehum coil) that is part of the refrigeration system. Dehumidification is done by utilizing an air pump. The pump injects fresh air into the chamber, displacing the humidified air. The pump cycles on/off based on the humidity setpoint. This results in dehumidification and humidity control without the use of a refrigeration system wet coil (dehum coil).

Refer to section Changing Humidity Setpoints to set the desired humidity setpoint.

LED Lighting (LGHT324, LGHT340, LGHT341, LGHT342)

White LED lights are standard on models 7350-25 and 7350-33, but are optional on models 7340-25, 7340-33. Colored LED lights are optional for both models.

Lights are controlled through the Light Cycle menu.



Model 7350

Operation of the Day/Night Setpoints

The system is used to simulate day and night testing.

This is a 24 hour cycle that can be programmed with separate temperature and humidity setpoints that correspond with the lights on (day) and lights off (night). The setpoints can also be made the same if continuous conditions are needed throughout the light cycles.

This option is broken down into 2 sections which work together as Day/Night simulation.

a) Controlling parameters and b) controlling lights.

a) Controlling parameters, Day/Night setpoint screen for parameters, daytime setpoint.




b) Controlling lights.

Each LED light bank has light bars mounted on each side of the chamber walls.




The Light Cycle screen has parameters that can be set up for the Day/Night simulation.

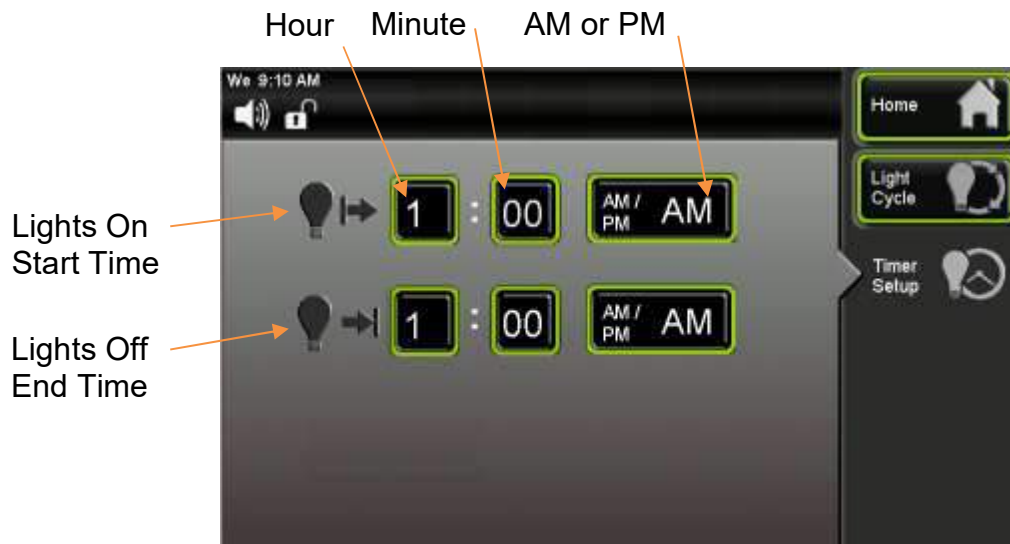

The Timer Setup screen allows users to setup lights “on” start time and lights “off” end time.


The Continuous Light button can also be used when the light cycle is setup for the light timer, this feature will disable any programmed Lights Start Time and Lights Stop Time.

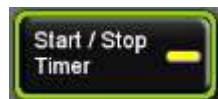

When the Start / Stop Timer (button) is active, the light timer will run the cycle based on the parameters that were setup. An icon appears in the status bar letting the user know that the Light Cycle is running (during this cycle the Continuous Light feature is disabled).



Continuous
Light

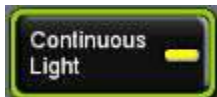

The Continuous Light (button) lets the light banks be on all of the time. When this feature is enabled, the features of the time based light cycle are disabled.

Setting the Light Cycle Timer




Lights
On/Off
Status

Light
Cycle
Menu


Screen shown with optional HUMD307,LED lights



To set the Light Cycle, press the (Light Cycle) button on the right side of the screen.

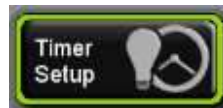

To set up the Light Timer press the (Timer Setup) button on the left side of the screen.

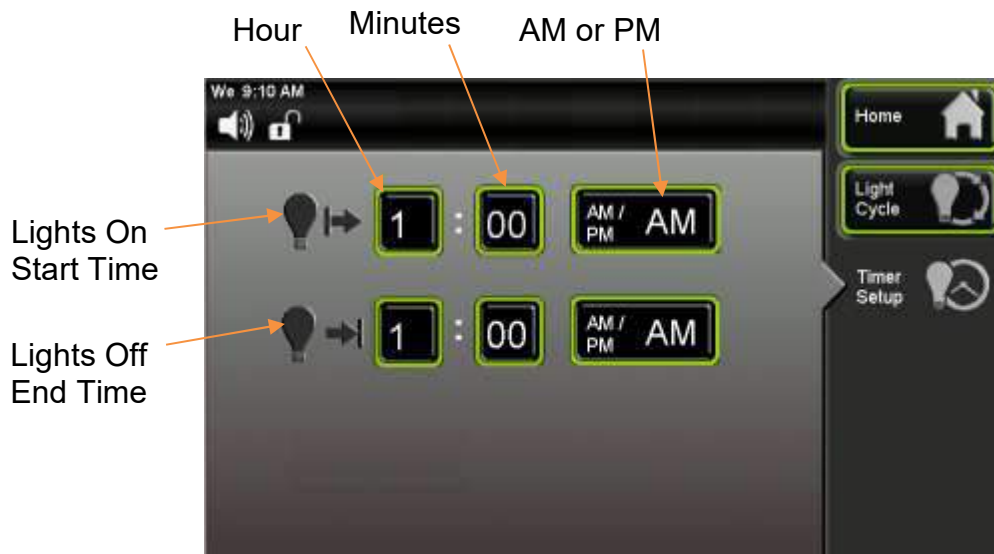

Press the Lights On Start Time hour button and the enter the hour that you want the lights to come on.

Then press the “minutes” button to set the minutes.



To setup AM/ PM, Press (AM /PM) button and the words for AM and PM will toggle back and forth.

Repeat process for setting up Lights Off End Time.



Go back to the previous screen by pressing (Light Cycle) button.



To activate the Light Cycle Timer press the Start / Stop Timer button.


Status of the Light Cycle


The light cycle is programmed to automatically change lights, temperature and humidity (optional) over a 24 hour period. The Light Cycle Info screen is intended to provide a comprehensive visual illustration of the Light Cycle settings and current conditions.

When the Light Cycle is running, the status of the cycle time can be displayed by pressing the Light Cycle icon that is in the Status Bar.




The Light Cycle Info screen displays information regarding the light cycle, start time, end time, light banks etc. This icon is only displayed in the Status Bar when the Light Timer is active and during the “lights on” time.

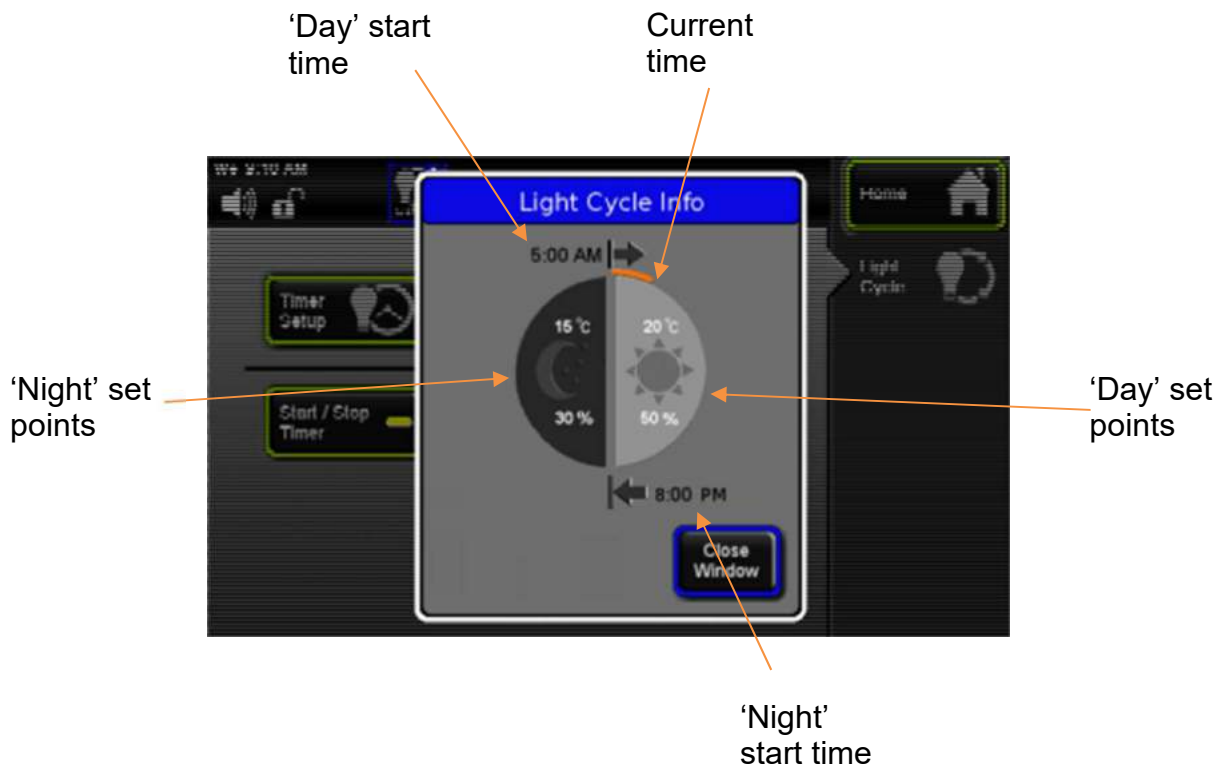

When Light Cycle is active the  Light Cycle Info icon appears in the Status Bar. Press this icon button and Light Cycle Info screen will appear displaying information about the Light Cycle and if it is near completion of the cycle.




To return to the home screen press the  Close Window button, or wait 15 seconds and the screen will return to the home screen.

Operation of Front Mounted 6" Recorders (RCDR316, RCDR317)

Built in 6" ink pen temperature and or humidity recorders can be purchased with CARON chambers. The recorders are shipped installed on the outer door of the chamber from the factory and require no installation.



Changing the chart paper:

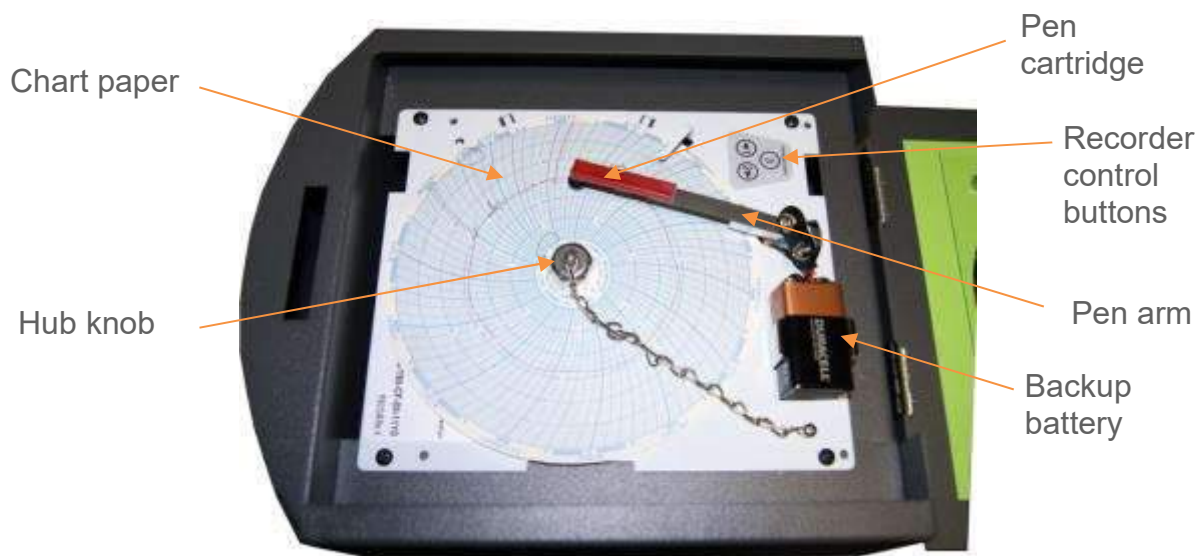

Press and hold the “change chart” button on the recorder (#3) for approximately one second until the pen begins to move to the left of the chart and then release the button. Wait until the pen has completely moved off of the chart. To remove the chart paper, unscrew (counter-clockwise) the chart “hub” knob at the center of the chart. Remove the old chart paper and position the new one so that the correct line coincides with the time line groove on the chart plate.

Re-attach the chart “hub” knob and fasten securely against the chart. Press and hold the “change chart” button (#3) again for approximately one second until the pen begins to move back onto the chart and then release the button. Check to make sure that the pen is marking on the chart paper. If it is not, then carefully adjust the pen arm to establish contact with the paper.


Chart recorder marking system:

This type of pen consists of a self-contained ink reservoir with a porous plastic stylus which is snapped around the outer edge of the metal pen arm. A pen cap is provided to extend the life of the ink pen during shipping or when the recording unit is not in service.

To remove the pen cap, gently lift the pen arm away from the chart paper. Remove the black plastic pen cap to expose the fiber tip of the ink pen and gently place the pen back onto the chart paper. Do not let the pen arm “snap” back onto the chart paper. This will flatten the fiber tip of the pen and will no longer give you a fine line marking on the chart paper. Place the pen cap in a safe place for future use. If the stylus does not touch the chart, adjustment can be made by slightly bending the metal pen arm in the center towards the chart paper. Do not use more pressure than is necessary to create a fine line marking on the chart paper. As the pen ink supply runs out, the pen color will become lighter. This indicates that the pen should be replaced.

Replacement of the Pen:

Recorders that are equipped with fiber tipped cartridge pens will have a cartridge that is color coded “red” to designate pen number one and an optional cartridge that is color coded “blue” to designate pen number two. The pen cartridge is securely fastened to the metal pen arm using a special “U” clip tab. For ease of replacement, it is suggested that the two screws that hold the pen arm be loosened and the pen cartridge and metal pen arm be removed as an assembly. Unsnap the plastic “U” clip tab of the pen cartridge from the metal pen arm, remove and discard the old pen cartridge. Replace the new cartridge by opening the hinge and snapping it securely around the metal pen arm. Refer to the image below:

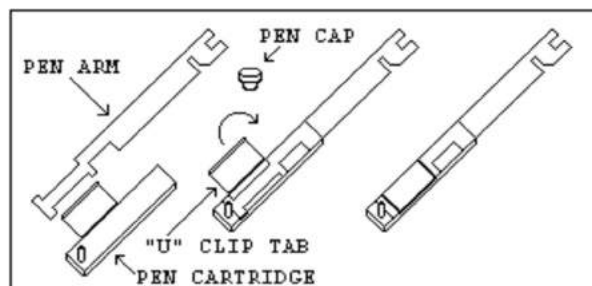

Pen Arm Calibration:

To check and/or adjust the recording pen(s) calibration to the outer most temperature graduation of the chart, press and hold the "change chart" button (#3) until the pen begins to move off of the chart. Once the pen(s) has moved off of the chart, again press and hold the "change chart" button (#3) until the pen begins to move back onto the chart. The pen should briefly stop at the outer most temperature graduation of the chart before continuing onto the chart to begin recording. If the pen does not stop exactly at this location on the chart, it can be adjusted or "calibrated" by using the left (#1) or right (#2) arrow buttons.

When the pen moves back onto the chart and briefly stops, you will have approximately five seconds in which to adjust the pen's position using the left and right arrow buttons of Figure 3.

On multiple pen recorders, each pen will move (one at-a-time) onto the chart briefly stopping at the outer most temperature graduation of the chart at which time the pen's position can be adjusted by using the left (#1) or right (#2) arrow buttons. When the time to adjust the position of the first pen has expired, the second pen will move onto the chart briefly stopping at the outer most temperature graduation of the chart at which time the second pen's position may be adjusted.

Each time the chart paper or fiber tip pen cartridge is changed, you should make sure that each pen stops at the outer most temperature graduation of the chart paper. Otherwise, this pen offset will cause the unit to record an incorrect temperature on the chart.

Recorder Calibration:

If a calibration adjustment is required for a single pen recorder, use the left (#1) and right (#2) arrow push buttons on the recorder to calibrate (or move) the pen's position on the chart to correspond to the temperature of the solution. The arrow buttons must be held for approximately five seconds before the pen will begin to move.

For two pen recorders, you must first select the pen that you wish to calibrate. This is done by pressing the left (#1) arrow button to select the red pen or the right (#2) arrow button to select the blue pen. The arrow button must be held down until the green LED light goes out. After the green LED light goes out, follow the instructions in step #3 above.

Battery Backup:

The green LED light remains a constant green color indicating that both the battery and the main power to the unit are good. Refer to Figure 5 for the location of the green LED indicating light. If the AC power were to fail or the battery becomes weak, then the green LED light will begin "flashing" indicating that either you have lost the main power to the unit or it is time to replace the battery. Having a 9 volt DC battery back-up in place, will allow the recorder to continue to function normally for approximately 24 hours in the event of a power failure.

Operation of Front Mounted 10" Recorders (RCDR318, RCDR319)

Built in 10" thermal pen recorders can be purchased with CARON chambers. The recorders are shipped installed on the outer door of the chamber from the factory and require no further installation. Unlike ink pen recorders, the thermal recorders draw their own chart and control lines.

The 10" recorders have been setup at the factory in the following configuration: 7 Day / 24 Hour / Temperature 0-100°C / Humidity 0-100% (for dual input recorders). If this is not the ideal configuration for an application, the recorder may be reconfigured using the following process:

Configuring the recorder:

In order to configure the recorder, you will need to enter the set-up mode of the recorder. To enter the set-up mode of the recorder, press and hold the Change Chart button (#3) until the thermal pen arm begins to move off scale and then release the button. Similar to the 6" recorder.

Note: The green LED light will flash fast while the thermal pen arm is moving off scale.

Wait until the thermal pen arm has moved completely off scale and stops (the green LED light will stop flashing and will be steady On). Unscrew (counter clockwise) the chart "hub" knob at the center of the chart and remove the recording chart paper. Gently lift the thermal pen arm just enough to be able to slide the paper out from beneath it. Remove the recording chart paper and place the Setup Chart onto the recorder. This chart contains the configuration categories of the recorder (Probe Input, Inner Chart Temperature, Outer Chart Temperature, Temperature Scale, Chart Rotation Speed, Input Filtering, Optional Relay Contacts and Date/Time for internal clock).

Next, press and hold either button #1 or #2 until the green LED light goes out and release the button. If this step is successfully completed, the pen arm will move to the outermost graduation ring of the Setup Chart. Use the Left (#1) or Right (#2) arrow buttons to adjust the center of the thermal pen to be on this outermost graduation ring.

Position the Setup Chart so that the tip of the thermal pen is in the center of the Start circle. Tighten the chart hub knob to secure the chart in place. Next, press and release the Change Chart button to begin. The chart will rotate to the first category (Input #1). Use the Left and Right arrow buttons to move the thermal pen arm to the desired option of each category. Press and release the Change Chart button to accept the selection and advance to the next category. You must press and release the Change Chart button when you have finished configuring the last category in order to save all of the changes that have been made to the recorder's configuration.

The thermal pen arm will move off of the chart allowing you to place the recording chart paper onto the recorder. Press and release the Change Chart button to begin recording.

Changing the Chart Paper:

Press and hold the Change Chart button (#3) for approximately one (1) second until the pen begins to move off scale and then release the button.

Note: The green LED light will flash fast while the thermal pen arm is moving off scale.

Wait until the thermal pen arm has moved completely off scale and stops (the green LED light will stop flashing and will be steady "On"). To remove the chart paper, unscrew (counter clockwise) the chart "hub" knob at the center of the chart. Gently lift the thermal pen arm just enough to be able to slide the paper out from beneath it. Remove the old recording chart paper and position a new one.

Re-attach the chart "hub" knob and screw securely (by hand) against the chart. Press and hold the Change Chart button (#3) again for approximately one (1) second and the thermal pen arm will move back onto the chart and begin recording.

Green Light LED Status:

The green LED light (located just below the three button membrane switch) is used to show the recorder's status:

- 1.) LED on steady (not flashing) and input(s) recording within chart range, indicates unit is recording normally.
- 2.) LED on steady (not flashing) and pen arm above outermost graduation and not moving, indicates recorder is in Change Chart mode. Press and release Change Chart button to return to normal recording mode.
- 3.) LED flashing rapidly and one or both inputs recording at outermost or innermost graduation indicates a sensor break. Check or replace sensor(s). If sensor(s) are ok, make sure process temperature is within configured range of recorder.
- 4.) LED flashing slowly (.8 seconds ON / .8 seconds OFF) indicates recorder is in Set-Up mode. Refer to section CONFIGURING THE RECORDER.
- 5.) LED is Off indicates that there is no power to the recorder. Check A/C power to the recorder.

Recorder Calibration:

If calibration is required for single input recorders, use the Left (#1) and Right (#2) arrow buttons on the recorder to calibrate the temperature being recorded on the chart to correspond to the temperature of the solution. The arrow buttons must be held for approximately eight (8) seconds before the pen begins to move.

If calibration is required for dual input recorders, you must first select the input that you wish to calibrate. This is done by pressing and holding the Left (#1) arrow button to select Input #1 or the Right (#2) arrow button to select Input #2. The arrow button must be held down until the green LED light turns off, after which follow the instructions in single input instructions above.

Maximizing Pen Life:

In order to maximize the amount of life expected out of the thermal pen tip, follow these simple rules:

- 1) Never let the thermal pen tip ride on the chart plate when the chart paper is not present. This will damage the protective coating of the heating element.
- 2) Never use chart paper that is creased or that has been folded.
- 3) Periodically clean the thermal pen tip with a cotton swap dipped in alcohol. Clean more often when operating the recorder in a dusty environment.
- 4) Always keep the door closed while the unit is recording.
- 5) Never lift the pen arm more than is necessary to remove and replace the chart paper. Excessive lifting may cause a decrease in the pen tip pressure and cause light printing.


Operation of H2O2 Sterilization cycle (STER305)

If the incubator is equipped with a STER305, then it is able to use the STER301 sterilization module (not included). STER305 includes an internal wiring connection for the sterilization module as well as door safety interlock. See separate STER301 User's Manual for details. Abide by all warnings.

Operation of Ramp/Soak Control (RAMP301)

A ramp and soak control system allows the user to store up to 50 steps spanning 5 profiles. A step consists of a change in setpoint (ramp), or maintain a setpoint for a fixed duration (soak). A step(s) can be repeated any number of times from 1 to 0 255 or infinite (repeat).

There are up to 4 different parameters that can be set up for the Ramp/Soak control: Temperature, Humidity, CO₂ and Lights (select models & accessories).

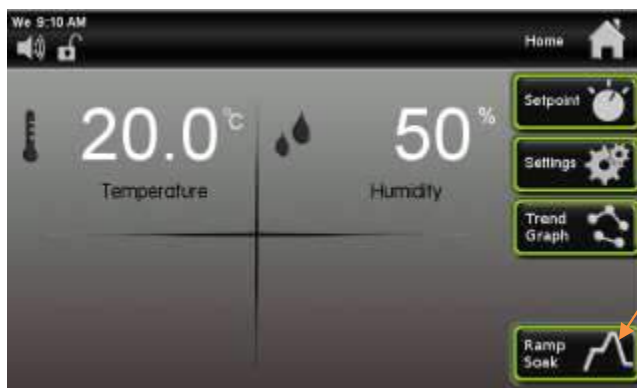

Ramp Soak
button



Press the (Ramp Soak) button on the Home screen.

Profile
button



Start / Stop
Ramp profile
button



Press the (Profile) button.

Profile Step buttons



Current Profile

Continue to Profile button



Press the (Step 1) button.

Ramp button


Unused button



Repeat button

Current Step

Soak button

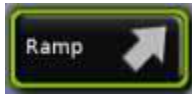

Press the (Ramp) button.

Ramp Rate parameters



Parameters that can change for profile

Step type

To set the ramp temperature value, press the (Temperature) button.



Enter the value of 30 and press (Enter).







Ramp rate allows a user to control the rate which parameters (temperature, humidity, CO₂, lights) change when transitioning to a new setpoint. This allows for a more gradual change than otherwise would occur.

Temperature tolerance range is 1°C when ramping temperature.

Humidity tolerance is 2% when ramping humidity.

There are (4) conditions to ramp to a new setpoint:

- 1) Ramp Rate = “0”, and “Wait for Process” button is not checked.
The chamber will work to reach the new setpoints as quickly as it can. The program will proceed to the next step immediately (before setpoints are reach).
- 2) Ramp Rate = “0” and “Wait for Process” button is checked.
The chamber will work to reach the new setpoints as quickly as it can. The program will wait until the chamber reaches the setpoint (tolerance) before proceeding to the next step.
- 3) Ramp Rate is a specified time, and “Wait for Process” button is not checked.
The chamber will work to maintain the ramp rate specified. The program will proceed to the next step when the time duration is complete, regardless if the setpoints have been reached or not.
- 4) Ramp Rate is a specified time, and “Wait for Process” button is checked.
The chamber will work to maintain the ramp rate specified. The program will proceed to the next step when the time duration is complete AND the chamber setpoints have been reached.

Using the time feature, the ramp cycle will run for the specified length of time then go the next profile. To set the ramp rate time, press the  (Hours), and or  (Mins) buttons.

Enter desired ramp time. (2 hours for example)



Wait for
Process



When (Wait for Process) button is selected, the parameter step will wait until the setpoints have been reached before proceeding to the next step .



Press the (Profile 1) button to go back to the Profile 1 screen to set up Step 2.

Step type
icon



Step type icon indicates what function will occur at that step.





Press (Step 2) button.





Press the  (Soak) button.




Press  (Hours) and or  (Mins) button(s) for desired soak time. This feature allows the unit to stay at the Ramp value for a specified amount of time before going to the next Step.

Enter values, example 2 hrs-10 mins.

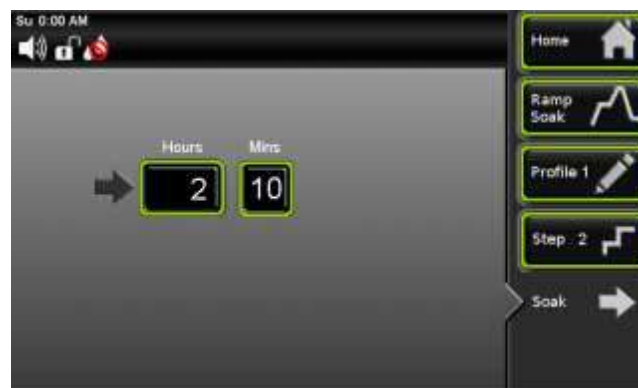



When finished entering values, press the  (Profile 1) button to go back to the Profile 1 screen to set up Step 3.




Press  (Step 3) button,



Choose the  (Repeat) button. To enter which step is to be repeated press the  and enter value on pop up key pad. There is a note that gives specific instructions which steps can be repeated. After entering the specified step, the number of repeats for that step can be specified. The # of repeats is a similar to a ‘loop’ that will allow a specified step to run continuously until the repeat cycle is complete, then it will move to the next step. The maximum number of repeats is 255.

What ***cannot*** be done is to create a repeat within a repeat or “nested” repeat. It is ok to have a repeat from step 7 to step 1 (black arrow), but not repeat step 9 to step 1 and include step 7 to step 1 (red arrow).

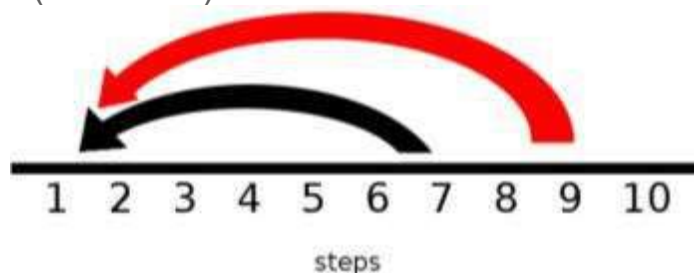



Enter values, example 2 for Step and 10 for # Repeats.




Once this is complete press the (Profile 1) button to return to the Profile 1 screen.



Press (Step 4) button.



Choose the Unused button. The Unused step is set up to be the “default” condition.

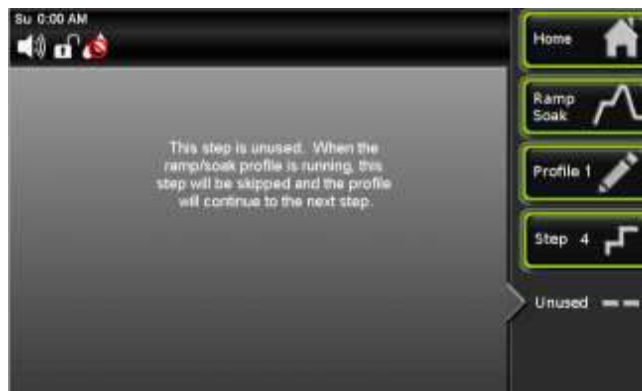

So when the Ramp, Soak or Repeat condition is not chosen for a step, the ramp profile program will skip over those unused steps.

When all of the Ramp profiles are entered, press the Start / Stop button and the ramp cycle program will run. The program can be turned off at any time during the process, but the Profile will start over from the beginning.

The other start / stop buttons for the other profiles will be greyed out indicating that only profile 1 has steps that can be run.

Ramp profile
running





The Ramp icon will appear in the status bar

Pressing this icon a pop up screen will appear with the profile steps that were chosen and what the current step the program is on. This is indicated by the orange bar.



When a ramp cycle has more than 10 steps, there is a button that will allow the steps to continue to the next profile. With the Profile screen open, at the bottom of the screen



Continue to
Profile
button



press the (Continue to profile) button. This will allow you to go to the next profile and set up the next number of steps. The numeric value shown will change based on the next available profile.



When finished creating of the required steps, press the (Ramp Soak) button to go back to Ramp Soak Profile screen.

Continue to
next Profile
icon



The black arrow indicates that after first profile 1 (1-10) is finished it will continue on to profile 2 (11-20).

This example screen is showing all 5 profiles are being utilized, up to 50 steps.



CALIBRATION

The temperature and optional humidity system can all be calibrated as necessary. CARON recommends an annual calibration check of each system. Before making a calibration adjustment, allow the cabinet to stabilize a minimum of 12 hours from a power off condition. If the unit has been in operation, allow a minimum of 3 hours of stable operation at all setpoints.


If you do not have the appropriate reference instruments to perform calibration, contact CARON's service department for on-site calibration at www.caronproducts.com Caron also provides validation services which ensure that the unit is functioning properly according to IQ, OQ and PQ protocols which satisfy FDA guidelines for qualification verification of equipment.




Be sure that all reference instruments are calibrated to an appropriate standard.

The Calibration Screen


To get to the calibration screen from the home page:

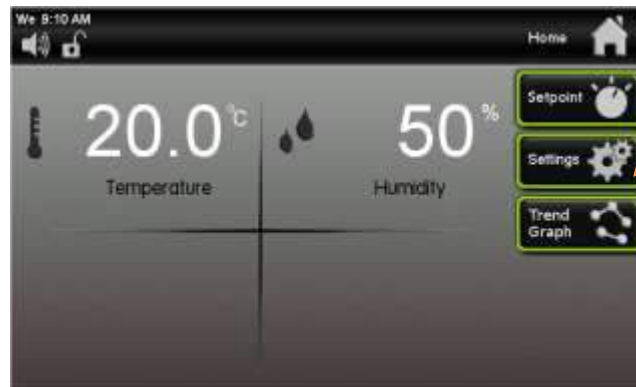


Settings
Button

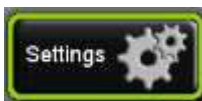

Press the (Settings) button.

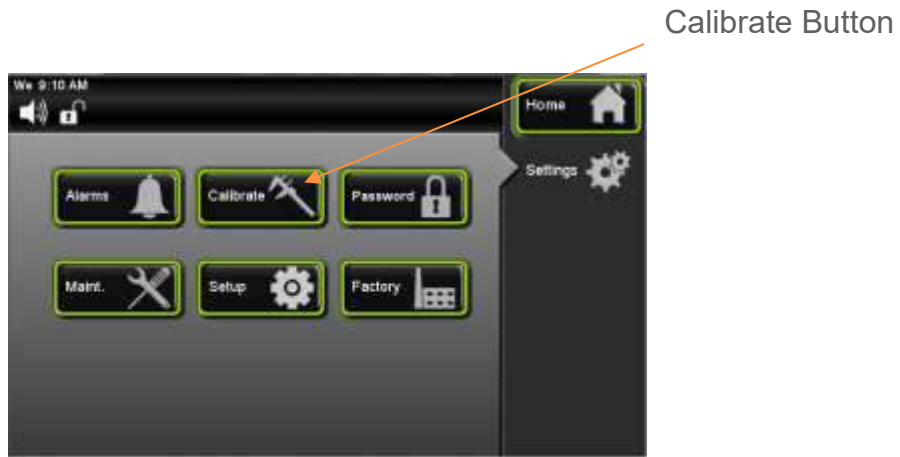

Press the  (Calibrate) button.

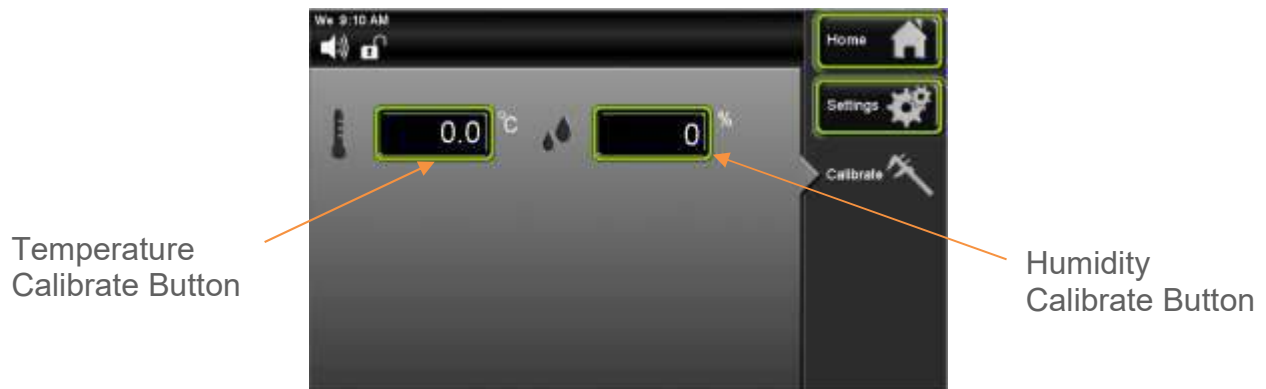


Calibration screen appears. If optional features such as Controlled Humidity are purchased, a calibration button will also appear for that option.


Calibrating Temperature

If temperature calibration is needed, the following steps can be taken:

Locate the reference instrument's temperature sensor in close proximity to the cabinet's geometric center. Be sure that the stabilization times described earlier have been satisfied prior to performing calibration.

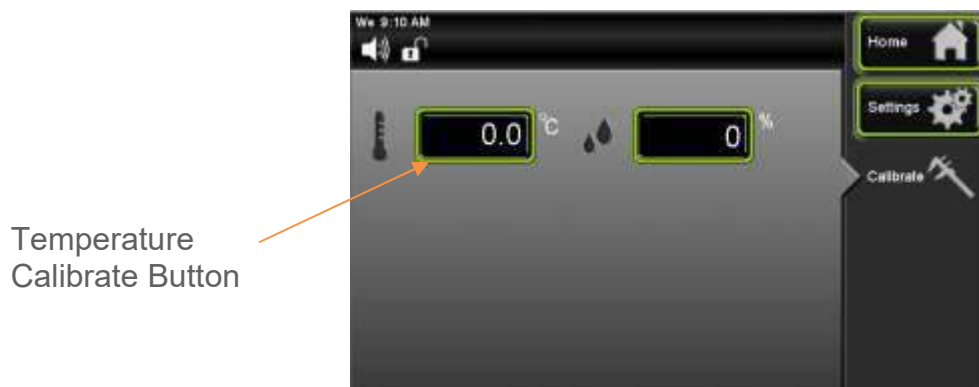

At the calibrate screen, press the  (Temperature Calibrate) button.

Enter the temperature offset by using the keypad and pressing  (Enter) when complete.

A positive value will move the temperature 'up' and a negative value 'down'. Press the 'home' button and verify the proper temperature is displayed.

Temperature calibration (example)

If the chamber temperature display reads 40.0°C and the calibrated independent sensor shows 40.3°C, set the temperature offset value to 0.3°C. If the calibrated independent sensor shows 39.6°C, then the entered offset should be negative. In this example the required offset to temperature would be -0.4°C.

Calibrating Humidity

If humidity calibration is needed, the following steps can be taken:

Locate the reference instrument's humidity sensor in close proximity to the cabinet's geometric center. Be sure that the stabilization times described earlier have been satisfied prior to performing this calibration.

A positive value will move the humidity 'up' and a negative value 'down'. Press the 'home' button and verify the proper humidity is displayed.

Humidity calibration (example)

If the chamber humidity display reads 80% and the calibrated independent sensor shows 83%, set the humidity offset value to 3.0%. If the calibrated independent sensor shows 74%, then the entered offset should be negative. In this example the required offset to humidity would be -6.0%.

Calibrating Optional Chart Recorders

For calibrating the optional front and side mounted chart recorders, refer to section (Optional Accessory Operation)

ALARMS

Alarm System Overview


The chamber control system is equipped with an alarm system that constantly monitors temperature, and humidity (on humidified models) to ensure the user is notified if the cabinet goes into an alarm condition. Notification occurs via an alarm pop-up window and a buzzer. Each alarm condition has been factory programmed to minimize nuisance alarms while maximizing warning time. There is a 2 hour time delay after start-up and setpoint changes. To avoid nuisance alarms after a routine door opening, an alarm condition must be present for 15 minutes* (45 minutes for humidity) before the operator is alerted. If the optional remote alarm contacts are present, in an alarm condition, the dry contacts will change state.

*Alarm delays are adjustable, see “Changing Alarm Setpoints and Delay” for details.


The following alarm messages may be displayed:

- Chamber temperature is higher than setpoint temperature
- Chamber temperature is lower than setpoint temperature
- Chamber humidity is higher than setpoint humidity (HUMD307 option)
- Chamber humidity is lower than setpoint humidity (HUMD307 option)
- Temperature sensor error

In the event an alarm occurs, the alarm indicator will appear on the status bar and an audible alarm pop-up window will automatically appear.

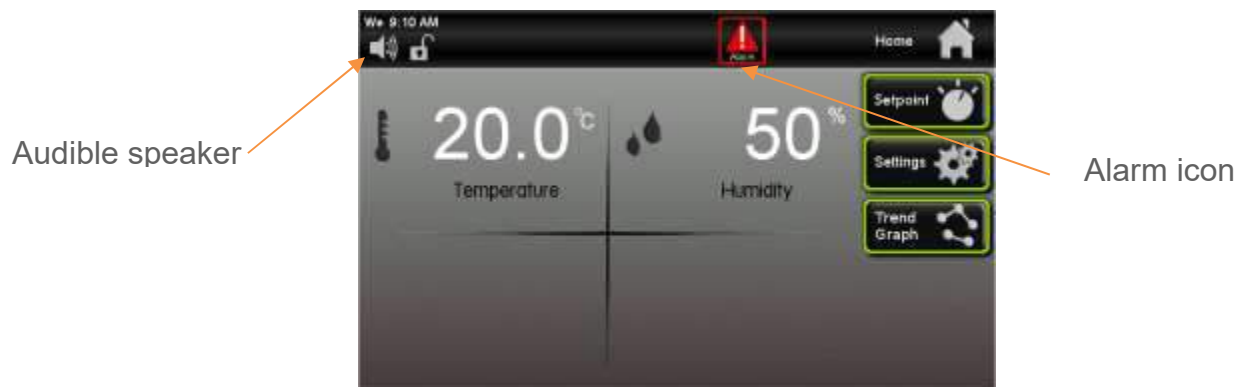


The flashing  (Alarm) icon will appear on the status bar.

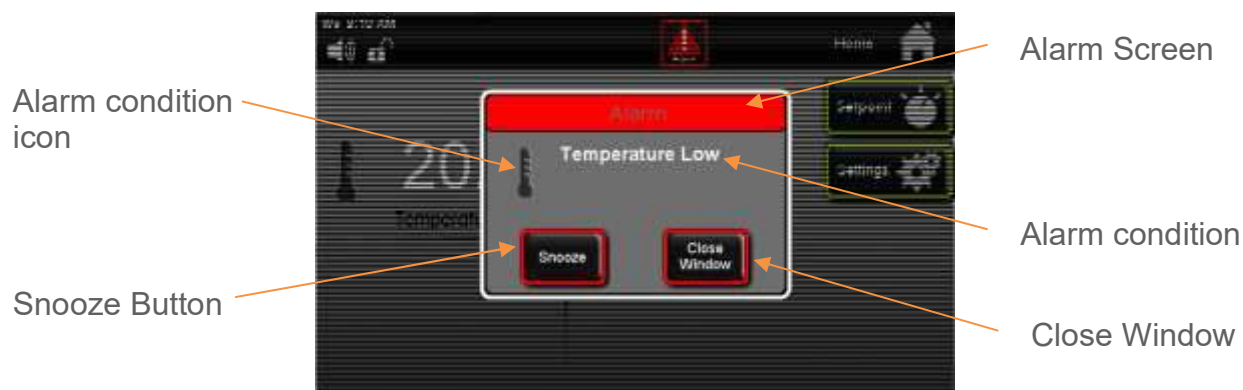

Audible Alarm Snooze Function:

When in an alarm condition, the Audible Alarm can be temporarily silenced to avoid being a nuisance to those nearby. The Audible Alarm will repeat after 1 hour has passed, if the condition has not been corrected. (The audible alarm will not sound if the alarm is muted, see Audible Alarm Mute)




Press the  (Snooze) button, the audible alarm is silenced for a period of 60 minutes.

When the alarm condition is corrected the alarm indicator and the audible alarm will automatically turn off (unless there is another alarm condition).




To check what the alarm condition is, press the  (Alarm) button on the status bar.



and the alarm window will be displayed. If the  (Snooze) button has already been pushed and 60 minutes have not passed the Snooze button will be “greyed” out.



If you press the  (Close Window) button, the Alarm Window will close, but the alarm will still be present as a flashing alarm icon on the status bar for the remainder of the 60 minutes time. It will not reset the 60 minutes alarm countdown time if the alarm condition is viewed on the pop up window.

After the 60 minutes time has passed for an alarm condition, the counter will reset itself to 60 minutes and repeat the countdown process again until the alarm has been resolved.

Audible Alarm Mute:


By factory default, when an alarm condition is present, the speaker will sound. This speaker can be muted in an 'on/off' fashion eliminating all audible sounds. (Muting the speaker will silence it until manually 'un-muted'. This is different than 'snooze' in the fact that snooze can only be enabled when an alarm condition is present and only lasts for 1 hour.) When the speaker is muted, the alarm icon continues to flash and the remote alarm contacts (optional) remain in the 'alarm' state.

To mute the speaker:





Press the  (Settings) button.



Press the  (Speaker Mute) button.




The Speaker Mute button toggles to the “on” position , and the speaker icon changes to  “Speaker Muted” icon.


Changing Alarm Setpoints and Delay

All alarm setpoints were pre-set at the factory to minimize nuisance alarms that could be created as a result of door openings. Alarm setpoints can be changed based on individual user requirements. Alarm values are deviations from the setpoint and are not actual setpoint values.

To change the alarm setpoints and delay:

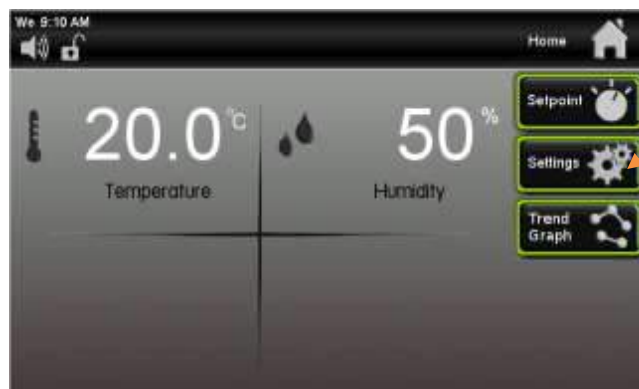

Settings
Button

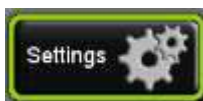

Press the (Settings) button.

Alarms
Button





Press the (Alarms) button.


Temperature
Alarm Button



Humidity
Alarm Button

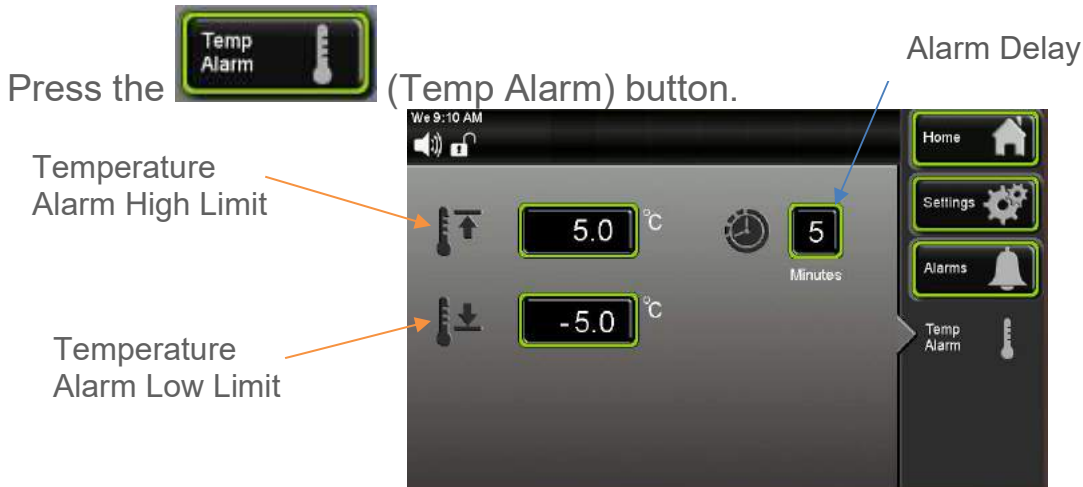




Once the alarm screen appears, press the  (Temp High Limit) button.

Keypad screen will appear. Enter the High Temp Alarm value; press  (Enter) when complete.

To change the Humidity Alarm, press the  (Alarms) button on the navigation menu to go back to the Alarms screen. Press the  (Humidity Alarm) button and repeat the same steps for humidity.

To change the alarm delay, press the alarm delay button  and enter the delay in minutes.

ALERTS

Alert System Overview


The chamber control system is equipped with an Alert system that constantly monitors features of the chamber and notify the user if the cabinet needs any type of service to ensure good running performance of the chamber. Alerts draw user attention to regular maintenance needs, and minimize the risk of a future alarm condition.


When an Alert notification occurs, contact www.caronproducts.com with the serial number of the chamber to order preventative maintenance kit(s).

Some of the Alert features are: Check the Atomizers (HUMD307 option), Replace the Air Filter, and Check Equipment Calibration is Due.

Notification occurs via an Alert icon on the status bar. When the Alert icon is pressed, a pop up window will display the alert condition(s). Each alert condition parameter is factory pre-set, no adjustment is necessary.



Press the  (Alert icon).

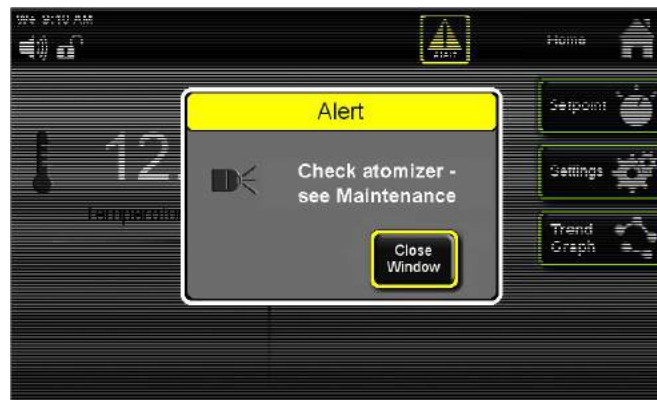



The Alert pop up window will appear displaying the alert message.




Press (Close Window) button to make the pop up window disappear.

Resetting Maintenance Alerts

Maintenance Menu Screen lets users check to see how much time is remaining on an item that may need routine service or calibration. This is very convenient to inform the user that a particular item will need to have service performed soon. After service has been completed, the item needs reset and the alert will disappear.

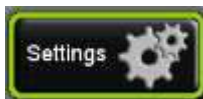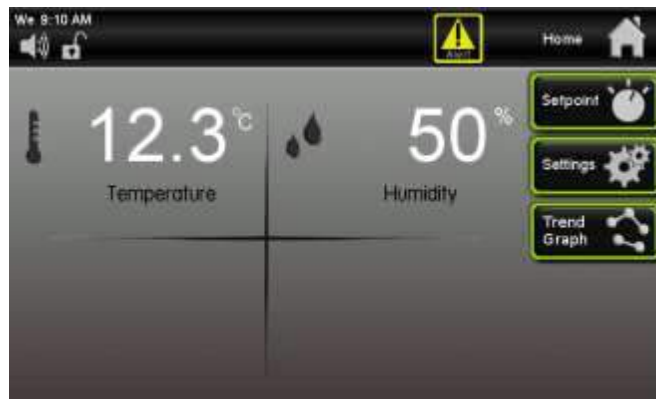

Press the (Settings) button.

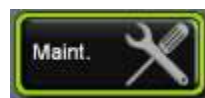

Once the Settings screen appears, press the (Maintenance) button.



Once a Maintenance item is displayed on the Alert screen, it will continue to be present as an icon in the Status Bar until the Maintenance item is corrected and the (Reset) button is pressed resetting the replacement time to “new” status.



Press the (Home) button to return to the main screen.


INFO

Info System Overview


The chamber control system is equipped with an Information system that constantly monitors the chamber and to notify the user when an automatic condition is occurring. This applies to conditions such as Defrost, DLOG301 or others that cannot be switched on and off by the user but is controlled automatically by the software of the control system. This notification cannot be disabled, it only lets the user know the chamber's current status.

Notification occurs via an Info pop-up icon on the status bar. When the Info icon is pressed a pop up window will display the Info condition(s).



The  (Info) icon will appear on the status bar.




Press the  (Close Window) button to return to the main screen.

ADVANCED FEATURES

Setting the Time & Day

The chamber has an internal real-time clock that keeps track of the day and time. It is set at the factory to Eastern Standard Time and may need to be adjusted for your time zone. To keep the clock accurate, it will need to be adjusted manually for daylight savings time changes. To set the day & time:

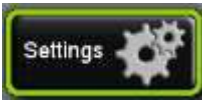

Press the (Settings) button.




Press the (Setup) button.




Press the  (Day / Time) button.




Press the  (Hour) button.

The Enter New Time in Hours window will appear. Enter the hour by using the keypad and pressing  (Enter) when complete.

Follow same procedure for setting up minutes.

To setup AM/ PM, Press  (AM /PM) button and the words for AM and PM will toggle back and forth.

To set the Day of the Week, press the  (Day of the Week) button. This button will scroll through the days of the week, press until the abbreviated letters correspond to the actual day of the week.

Press the  (Home) button to return to the main screen.

Door Heat

Some Caron units are equipped with an integral door heater to keep internal condensation off of the glass viewing area (not available on 7350). This is only for units with humidity control. In extreme ambient or running conditions, the factory default settings may need adjusted to maintain a clear viewing area. Increasing the door heater percentage will increase the amount of heat applied to prevent condensation. The door heater can run in either automatic or manual mode. Automatic is the preferred and default setting.



From the Setup screen press the  (Door Heater) button



In Automatic mode, the door heater works in conjunction with the internal heaters to maintain the temperature set point. In the event that the temperature is above set point, the door heater will automatically throttle back.

In this mode, the output percent is a scale factor of the overall heat output percentage. If condensation is present on the glass door under stable condition, then increase the door heater percentage value. The factory default value is 10%.



To change the output percent value Press the  (Door Heat Percent) button.

Enter the hour by using the keypad, press the enter button when complete.



In Manual mode, the door heater is a fixed output value regardless of temperature set point. This setting should only be used in units that have active cooling with temperature set points well above the low-end range.

Warning: in manual mode, the temperature set point may not be maintained.

Humidity Control (models with controlled humidity)

This feature allows the Humidity to be controlled by a setpoint determined by the user, which will be displayed on the home screen. If this feature is disabled, the humidity value on the home screen will be in a “read only” condition. An icon in the status bar will indicate whether the Humidity Control is enabled or disabled



From the Setup screen press the (Humidity Control) button


Humidity Control disabled icon



Humidity Control disabled indicator

When the humidity control is disabled the toggle button indicator is off and the humidity control icon in the status bar has a red circle around it.

Humidity Control enabled icon

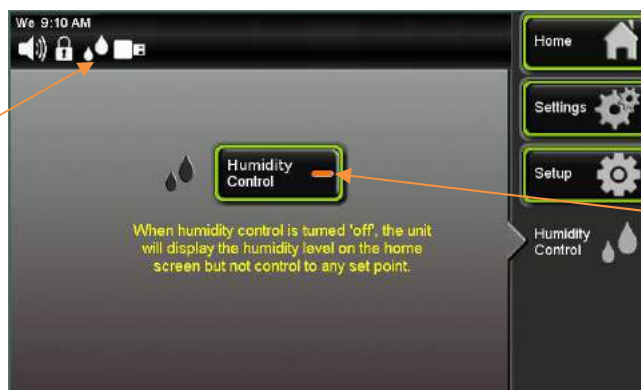


Humidity Control enabled indicator

When the humidity control is enabled the toggle button indicator is on and the humidity control icon in the status bar has the red circle removed. Once the selection has been made the status of the Humidity control will be visible on the home screen.

Locking the Controls

To prevent unauthorized and accidental changes, the touchscreen can be locked-out. The passcode is required to lock-out the controls and the same passcode is used to unlock it. The factory default passcode is '1234'. This passcode can be changed by the user to create a unique 4-digit passcode. There is also a feature that will let you change the passcode from the factory default to a user defined passcode. The factory default for the screen lock is "unlocked"

To lock the touchscreen,

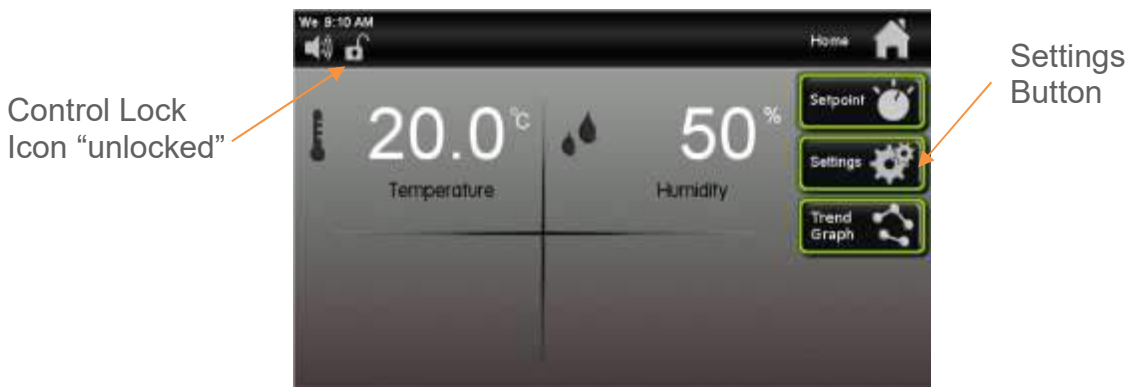

Press the  (Settings) button.



Press the  (Password) button.



Press the (Lock Keypad) button.




The Enter the Current Passcode keypad screen will appear.




Enter digits “1 2 3 4”; press (Enter) when complete.

The screen will change back to the Home Screen and the Control Lock icon will change to the “locked” position.

Control Lock
Icon "locked"

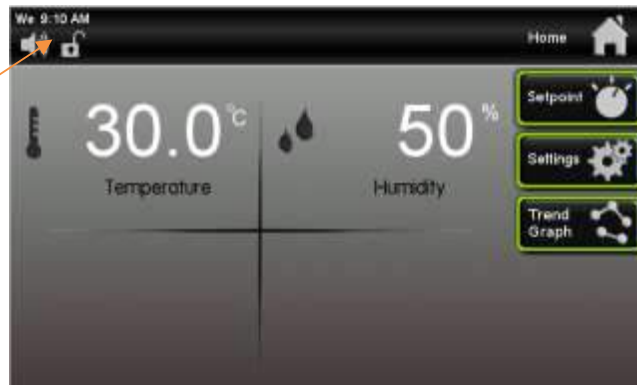

When any button is pressed on the home screen the following pop-up window will

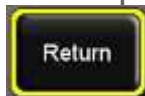

appear. If the button is pressed, the screen will change back to the Home Screen.




To unlock the touchscreen,




From the previous Alert "Keypad is Locked" pop up screen, press the (Unlock) button. The Enter New Passcode window will pop up.



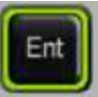

Enter the digits “1 2 3 4”; press  (Enter) when complete. The Control Lock Icon will change back to the “unlocked” position.


Control Lock
Icon “unlocked”

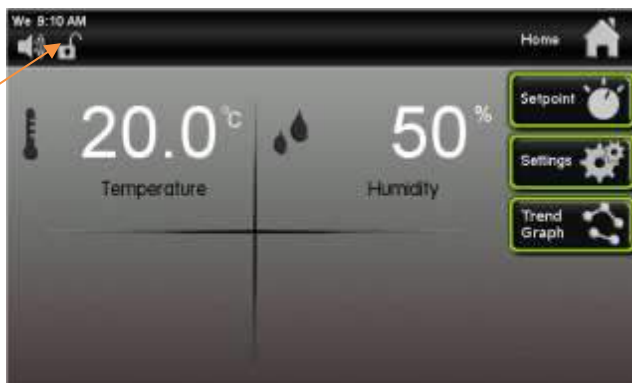

Changing the passcode

To prevent unauthorized and accidental changes being made to the chamber, the touchscreen can be locked-out. The passcode is required to lock-out the controls and the same passcode is used to unlock it. The factory default passcode is '1234'. This passcode can be changed by the user to create a unique 4-digit passcode. The current passcode is required to change the passcode.

To lock the touchscreen,

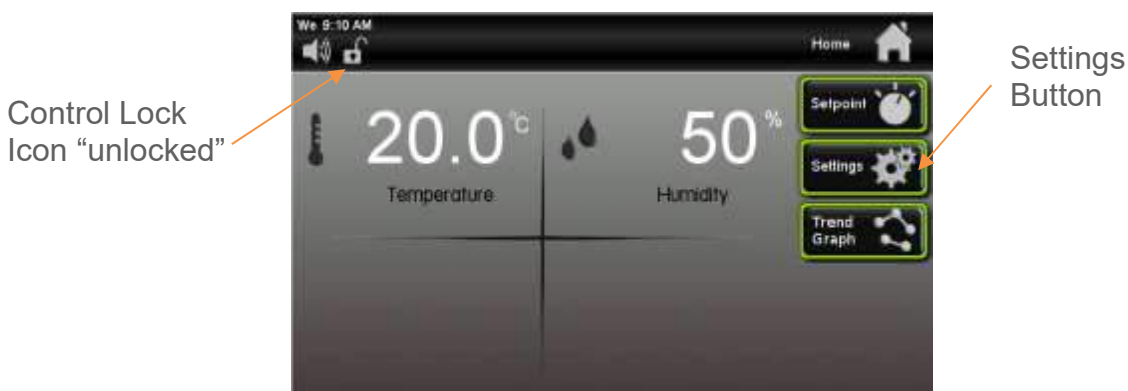


Press the  (Settings) button.



Press the  (Password) button.



Press the  (Change Passcode) button.



The Enter Current Passcode keypad screen will appear.



Enter digits “1 2 3 4”; press  (Enter) when complete.



The Enter New Passcode keypad screen will appear.

Enter any new four-digit passcode (example: “2 5 8 0”). Then press  (Enter) when complete.




The Lockout screen will tell you that the Passcode has been changed to a new value. *This is only time that the Passcode will be displayed on the Lockout screen.*

Factory menu & troubleshooting

The chamber control system is equipped with advanced diagnostics features which allow the user to manually turn 'on' & 'off' each electronically controlled system. The factory menu can be used to

- View the current chamber configuration
- See the percent output of the control system
- Manually and individually toggle any output

To access the Factory Menu,

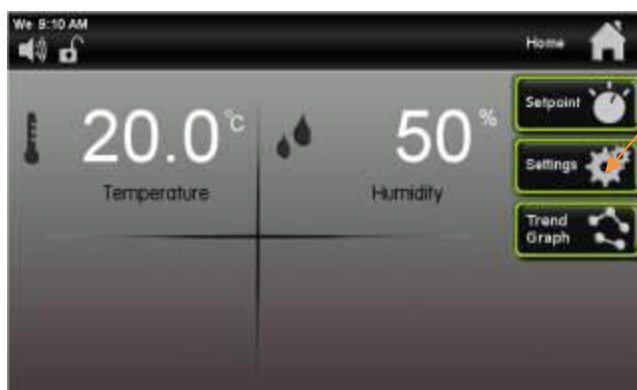

Settings
Button



Press the (Settings) button.




Factory
Button



Press the (Factory) button.



From the factory menu, four items can be selected. Press the (Model Details) button to view the chamber's configuration

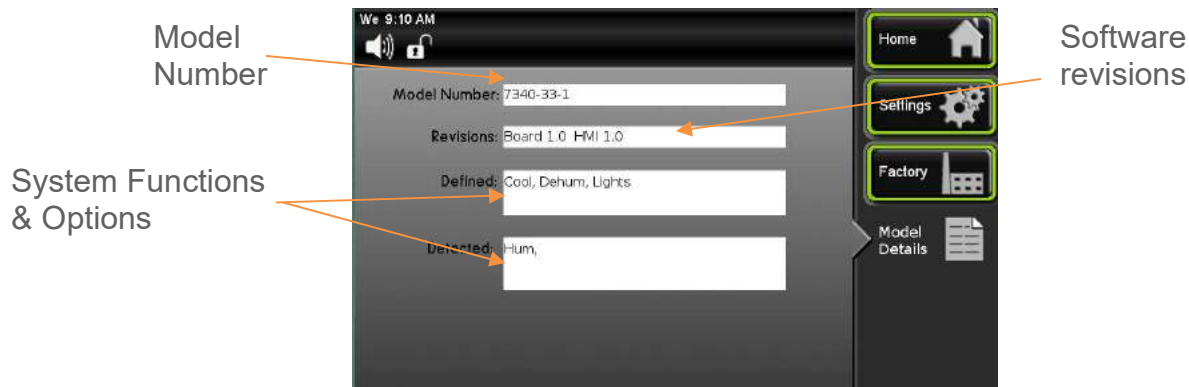

From the Factory screen, press the (Output Percent) button to view the current percent output level of each control parameter.




Output
1 & 2
Buttons



Navigating to the Output 1 or Output 2 screens in the factory page will temporarily halt chamber control & functionality.

To individually and manual control each output variable, from the factory screen press



the  (Output 1) button. Note: Based on the chamber model number and options, not all functions will be present.



Each item can be turned on to check the condition of that device or parameter to aid in diagnosing a problem.



Press the (Output 2) button for other parameter buttons.



Chamber control & functionality is restored as soon as the screen is exited (Home, Settings, or Factory buttons) .When finished with diagnosis in Output 1 or Output 2



screen, press the (Factory) button to return to that screen. Once you go back to the Factory screen all parameters that were selected in Output 1 or Output 2 screens will reset to the “off” position.



Press the (Home) button to return to the main screen.

PREVENTATIVE MAINTENANCE

The CARON chamber has been robustly designed to minimize performance problems. However, regular maintenance is very important for continuous trouble free operation.

As a general rule, CARON recommends an annual calibration check of the temperature and humidity systems. CARON offers a full range of on-site calibration and validation services. We also offer preventative maintenance contracts on our equipment. Contact our Service department for details at 740-373-6809 or visit us on the web at www.caronproducts.com.

Recommended Daily Maintenance Checks

- Check the Temperature and Humidity displays versus setpoints.
- Check for and correct any alarm condition.

Recommended Monthly Maintenance Checks

- Check to ensure the drain in the bottom of the unit is draining properly.
- Check front air intake filter. If the filter is dirty replace it with Caron Preventative Maintenance PM Kit. Washing the filter will result in poor performance.
-

Recommended Annual Maintenance Checks

- Replace humidity atomizer nozzle (HUMD311-1,-4 option) (see Replacement Parts section).
- Disinfect all interior surfaces with a general purpose laboratory cleaning agent.
- Perform a complete calibration of the temperature and humidity systems.
- A full validation is recommended for GMP facilities each time a unit is installed, moved or undergoes significant repair. Contact CARON's service department to schedule on-site validation.

Here is a list of PM Kits that are available for models and accessories covered in this manual.

| Model | PM Kit |
|--------------|---------------|
| 7340-25 | PM-7340-25 |
| 7340-33 | PM-7340-33 |
| 7350-25 | PM-7350-25 |
| 7350-33 | PM-7350-33 |

| Accessory | PM Kit |
|------------------|-------------------------|
| FLTR301 | PM-FLTR301 |
| HUMD311 | PM-HUMD311 |
| LGHT602 | ² PM-LGHT602 |
| RCDR316 | PM-RCDR316 |
| RCDR317 | ¹ PM-RCDR317 |
| RCDR318 | PM-RCDR318 |
| RCDR319 | ¹ PM-RCDR319 |
| RCDR320 | PM-RCDR320 |
| RCDR321 | ¹ PM-RCDR321 |
| RCDR322 | PM-RCDR322 |
| RCDR323 | ¹ PM-RCDR323 |
| STER301 | PM-STER301 |
| | PM-STER304 |

¹only if HUMD311 is installed

²only if LGHT602 is installed

SPECIFICATIONS

| MODEL | 7340-25 | 7350-25 | 7340-33 | 7350-33 |
|------------------------------|--|------------------|--|------------------|
| Temperature Range | 10°C to 40°C | | | |
| Temperature Control | ± 0.1°C | | | |
| Temperature Uniformity @20°C | ± 0.3°C | | | |
| Temperature Sensor | 3-wire RTD | | | |
| Lights | - | LED White | - | LED White |
| Evaporator coil | Electrostatic Coated Coil | Standard Coil | Electrostatic Coated Coil | Standard Coil |
| Fresh Air Ports | - | Standard | - | Standard |
| Interior Dimensions | 32" W x 27" D x 53" H (81cm x 69cm x 134cm) | | 32" W x 27" D x 66" H (81cm x 69cm x 168cm) | |
| Interior Construction | Type 304, 2B Finish, Solid Stainless Steel | | | |
| Exterior Dimensions | 36" W x 33" D x 77" H (90cm x 85cm x 196cm) | | 36" W x 33" D x 90" H (90cm x 85cm x 229cm) | |
| Exterior Construction | Cold Rolled Steel, Powder Coated | | | |
| Work Space | 25 Cu. Ft. (708 Liters) | | 33 Cu. Ft. (934 Liters) | |
| # of Shelves | 4 | | 5 | |
| Shelf Construction | Type 304, Perforated Stainless Steel, Electro polished | | | |
| Shelf Dimensions | 29" W x 24" D (74cm x 62cm) | | | |
| Door | Glass | Light Tight Door | Glass | Light Tight Door |

Specifications are subject to change without notice.

Environmental Conditions: Temperature 15°C to 25°C, Humidity non-condensing

*See graph for details


| Model | 7340-25,7350-25 | | | 7340-33,7350-33 | | |
|-----------------|--------------------|--------------------|-------------------------|--------------------|--------------------|-------------------------|
| | -1 | -2 | -3 | -1 | -2 | -3 |
| Shipping Weight | 575 lbs (261 kg.) | | 875 lbs. (397 kg.)** | 650 lbs. (295 kg.) | | 950 lbs. (431 kg.)** |
| Electrical | 115V, 60Hz, 16A | 230V, 60Hz, 10A | 230V, 50Hz, 8A | 115V, 60Hz, 16A | 230V, 60Hz, 10A | 230V, 50Hz 8A |

**Includes export shipping crate


*7340/7350-25 Series units have forced internal air flow of 350 cfm (9,900 LPM)

*7340/7350-33 Series units have forced internal air flow of 450 cfm (13,000 LPM)

ELECTRICAL SCHEMATICS

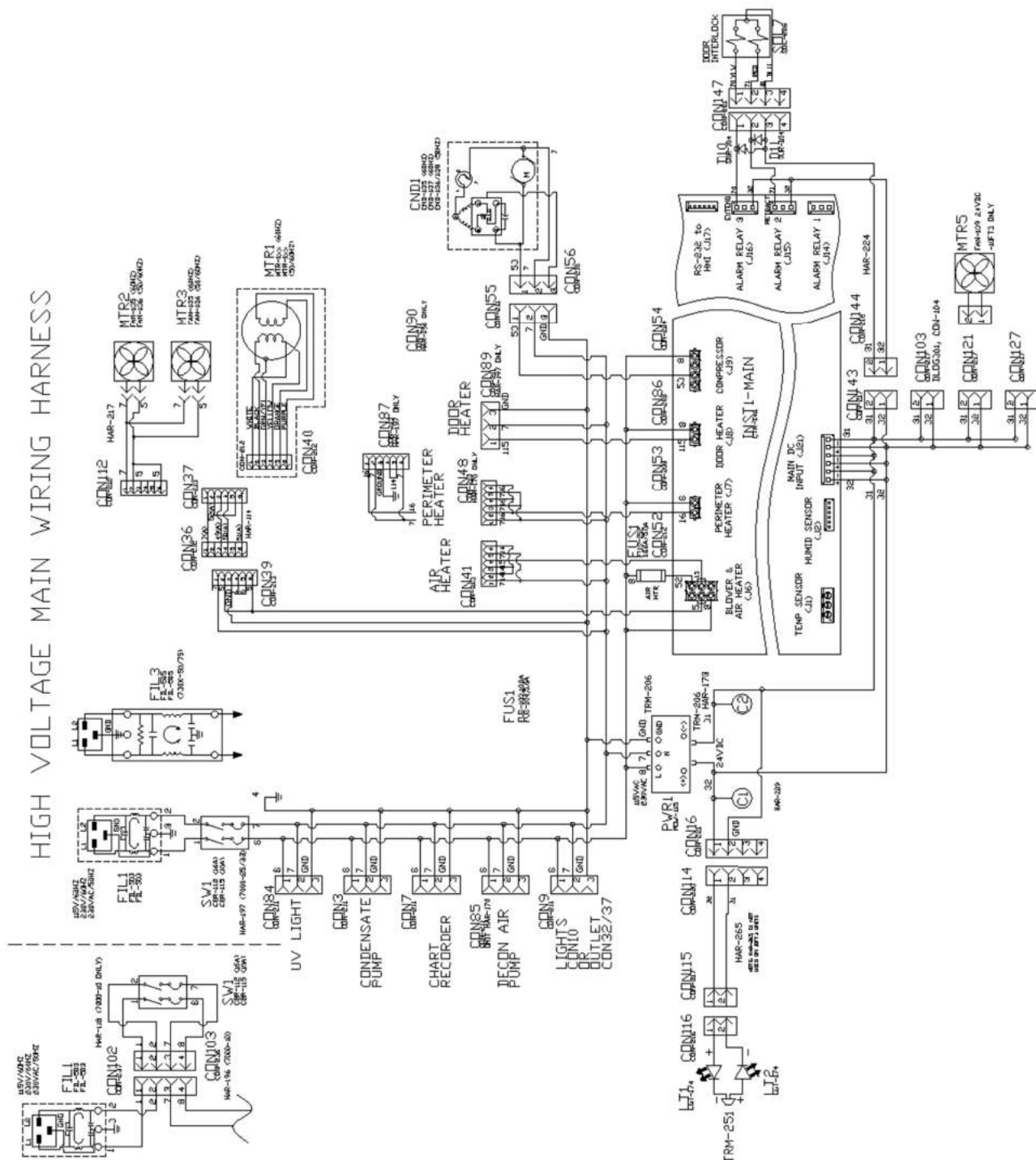

MAIN CONTROLLER BOARD (CTR-140)

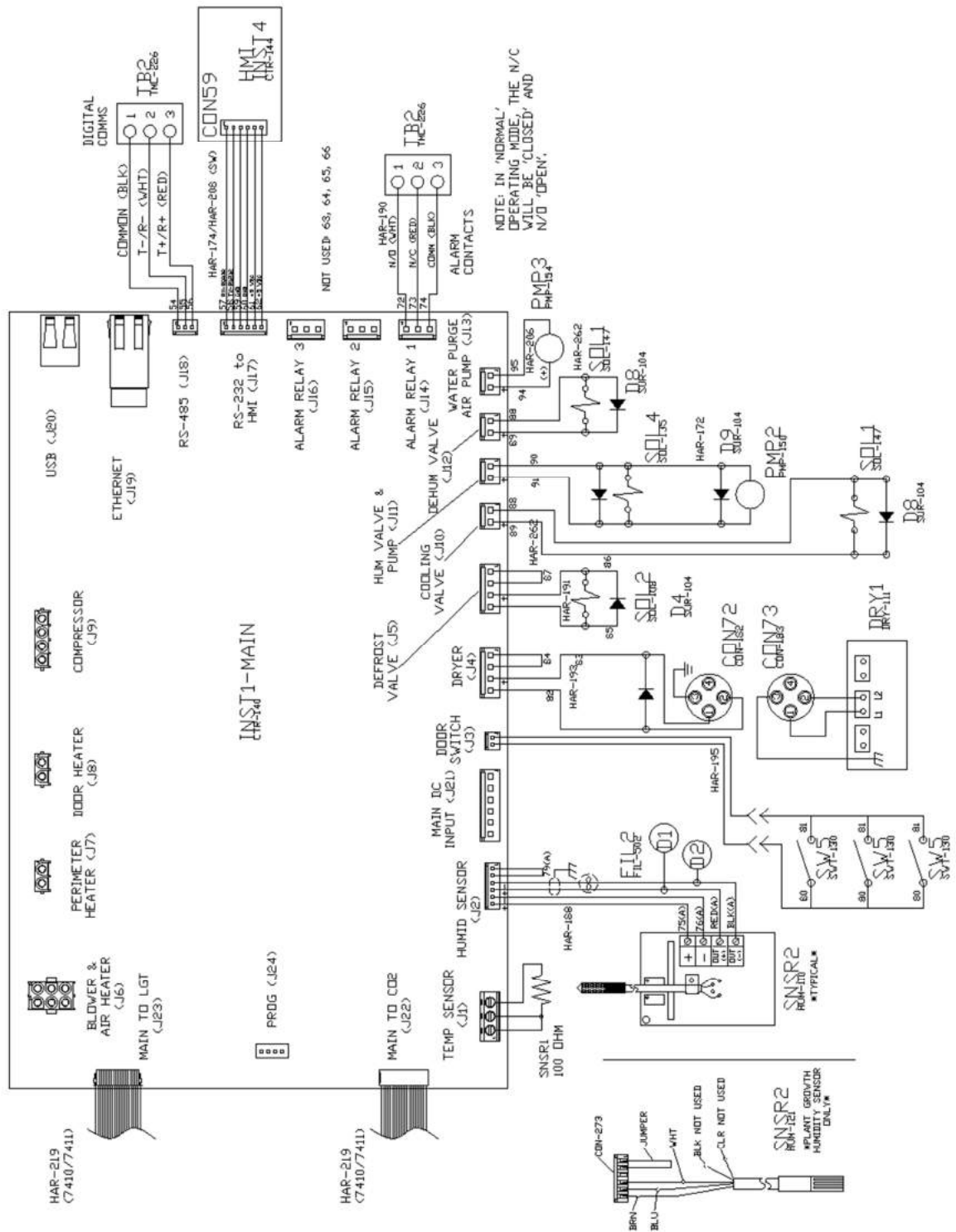

The diagram illustrates the wiring for the TORADIX IRIS system, showing connections between the main unit and various peripherals.


Main Unit (TORADIX IRIS):

- CON106:** Connector for the CTR-146, showing pins 31 and 32.
- MEM-104:** Memory module.
- POWER 24VDC:** Power input.
- CON107:** Connector for the HAR-214.
- USB:** USB port.
- ETHERNET (J19):** Ethernet port.

Peripherals and Connections:


- INST1-MAIN (CTR-140):** Main unit with connectors CON105, RS-485 (J18), RS-232 to HMI (J17), and ALARM RELAY 3 (J16).
- HAR-213:** Cable connecting CON105 to CON106.
- HAR-214:** Cable connecting CON107 to CON108.
- CON108:** Connector for the HAR-214.
- CON109:** Connector for the HAR-213, showing pins 54, 55, and 56.
- CTR-146:** Cable connecting CON109 to CON106.
- CON105:** Connector for the INST1-MAIN, showing pins 54, 55, and 56.
- RS-485 (J18):** RS-485 port on the INST1-MAIN.
- RS-232 to HMI (J17):** RS-232 port on the INST1-MAIN.
- ALARM RELAY 3 (J16):** Alarm relay on the INST1-MAIN.
- MAIN DC INPUT (J21):** Main DC input on the INST1-MAIN.
- DOOR SWITCH (J3):** Door switch on the INST1-MAIN.
- DRYER (J4):** Dryer on the INST1-MAIN.
- HAR-173:** Cable connecting CON103 to CON104.
- CON103:** Connector for the HAR-173, showing pins 31 and 32.
- CON104:** Connector for the HAR-173, showing pins 31 and 32.
- CON-217:** Connector for the HAR-173.
- CON-216:** Connector for the HAR-173.
- HAR-215:** Cable connecting CON106 to the USB Panel Mount.
- USB Panel Mount:** USB panel mount.
- MEM-103:** Memory module.
- HAR-216 TBD:** Cable connecting CON107 to the USB Panel Mount.

LIGHT CONTROLLER BOARD (CTR-142)

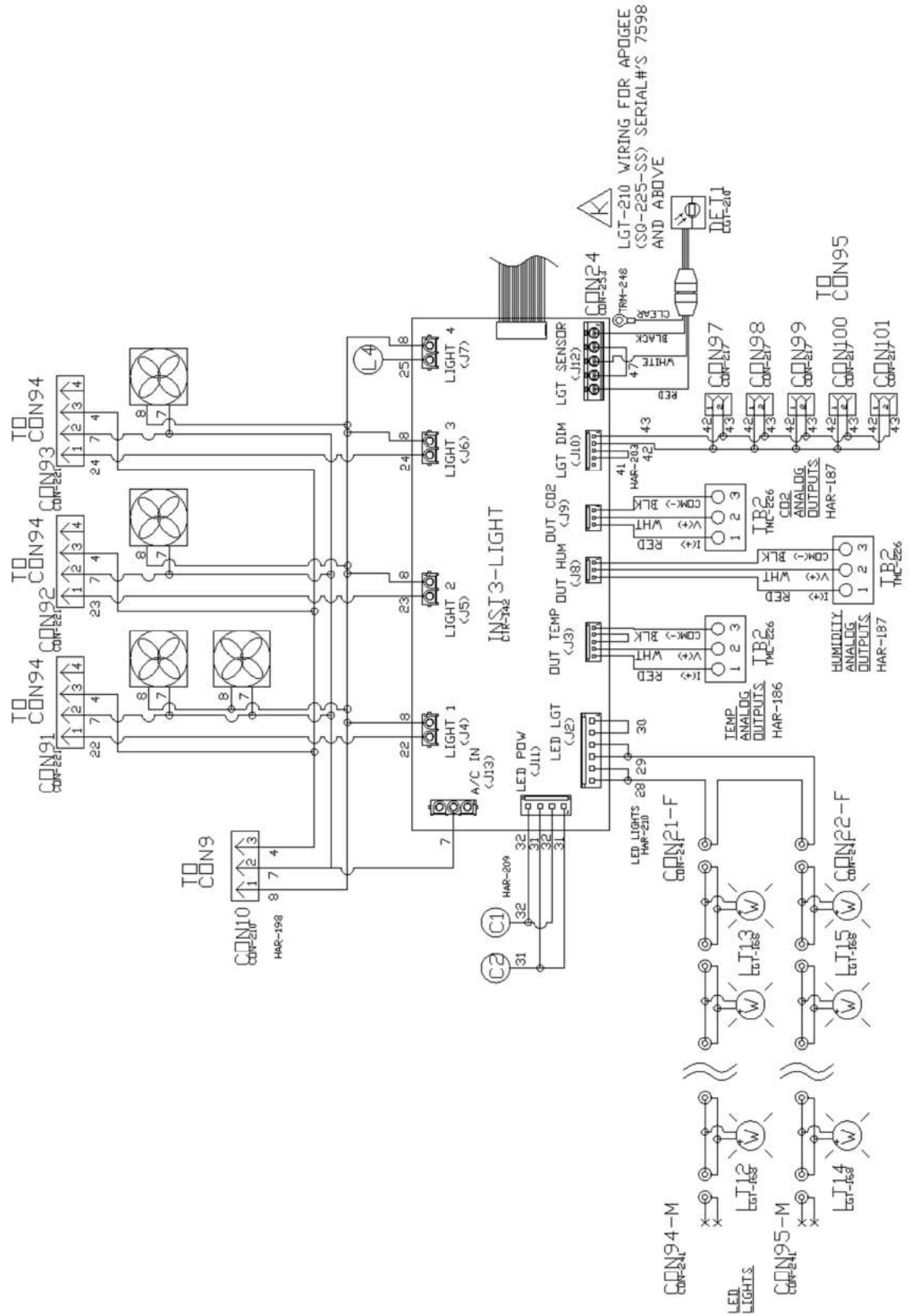


AIR HEATERS




7300-75-2/3, 7301-75-2/3,
7302-75-2/3, 7303-75-2/3
NO HAR-114 ADAPTER HARNESS
EVEN THOUGH IS 230V & 2X
HEATERS

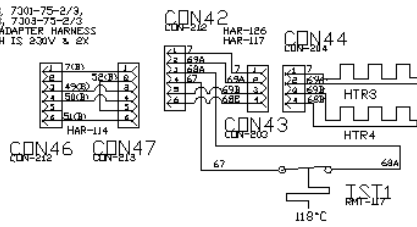


7304-75 FT³/3 PLANT GROWTH (1 HEATER)

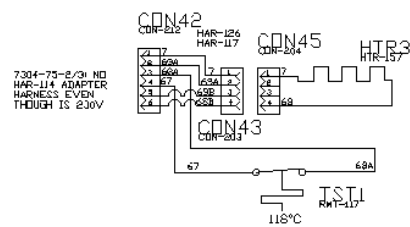


DOOR HEATER

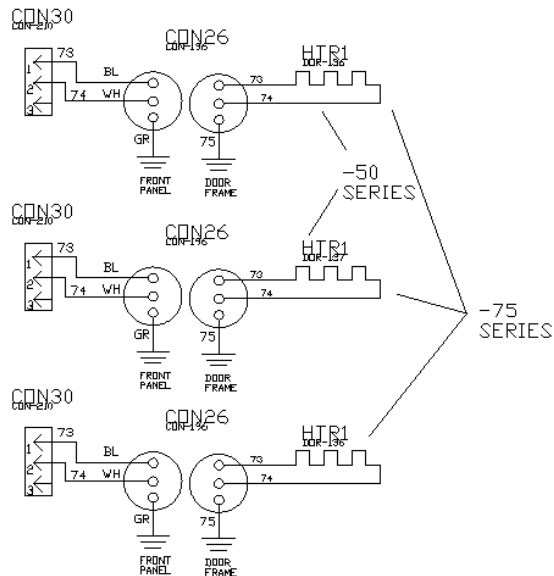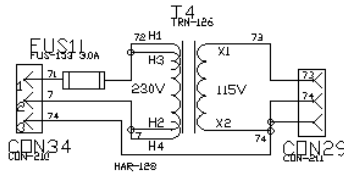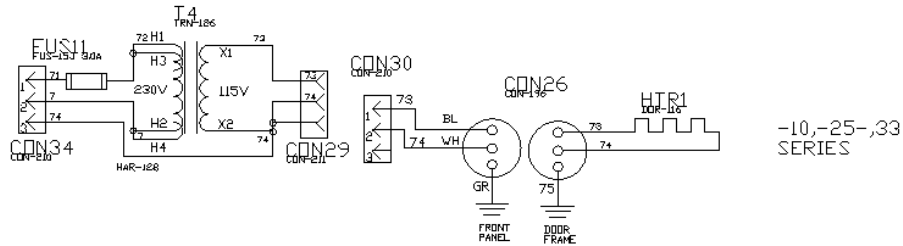



CO2 CONTROLLER BOARD (CTR-141)

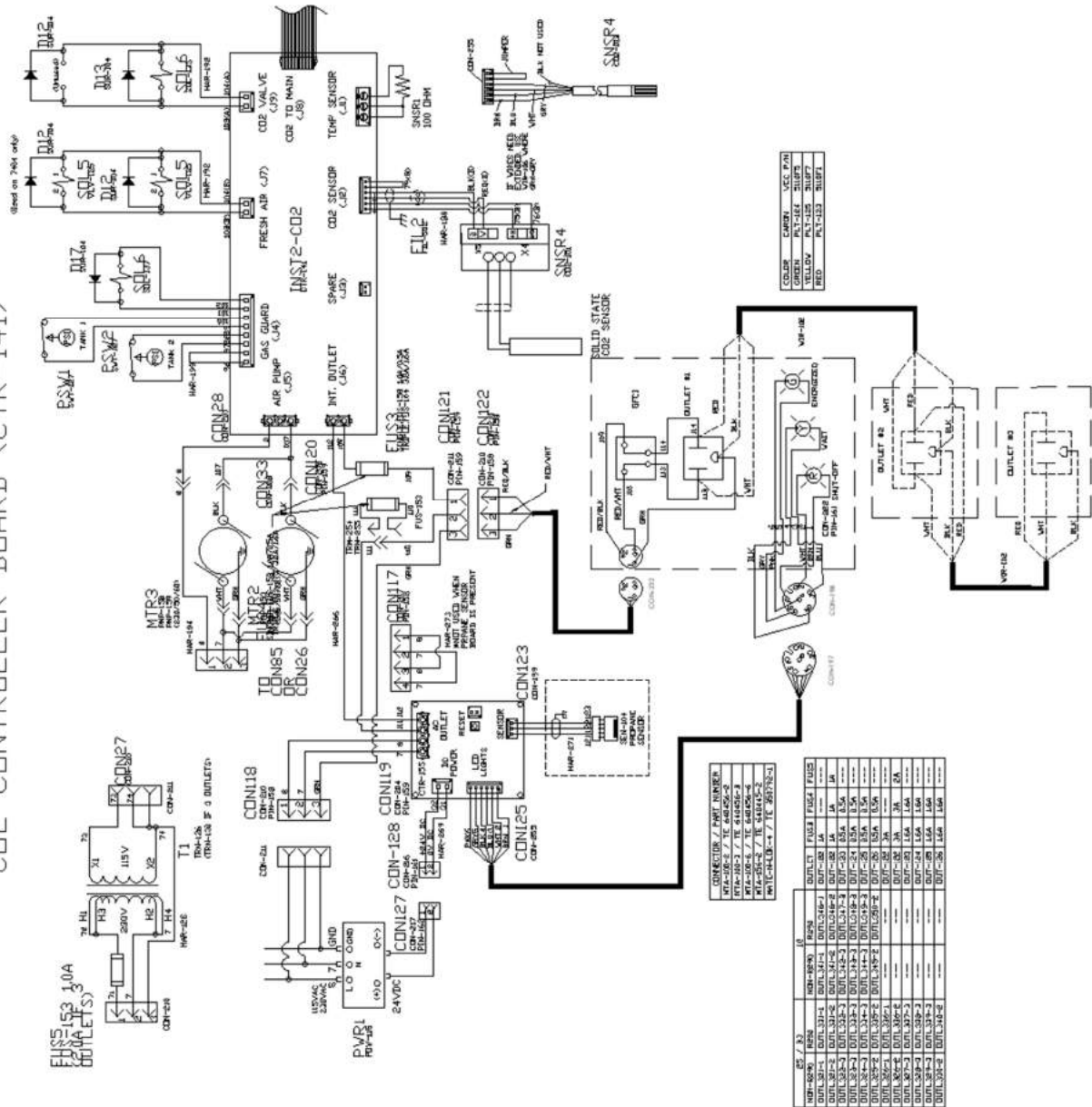


PERIMETER HEATERS

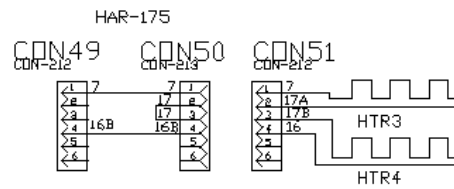


75 FT3 ONLY

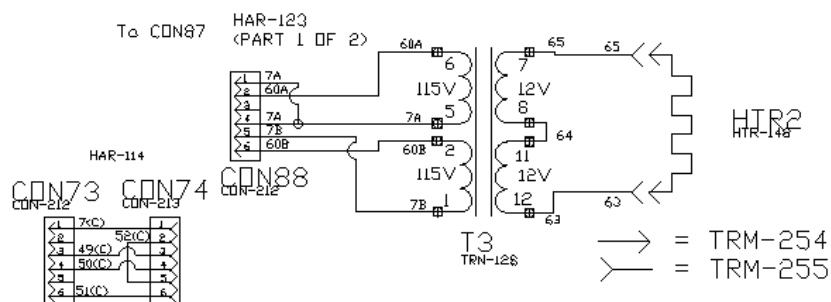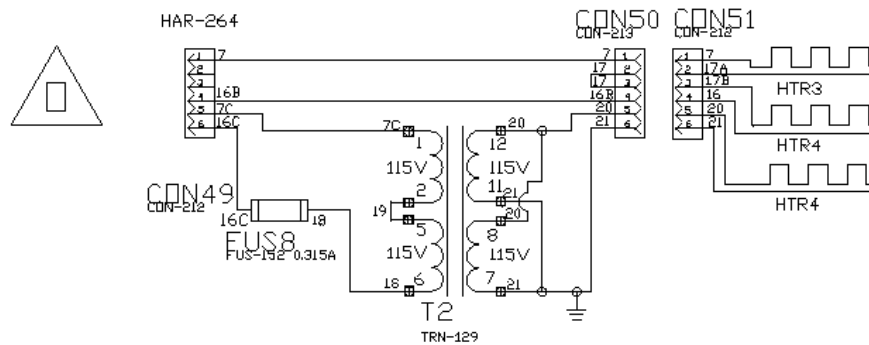


UV LIGHT

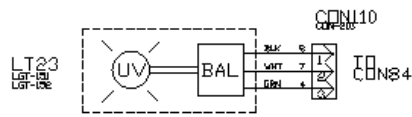

CONDENSATE PUMP

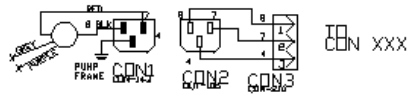


FRONT MOUNT RECORDER

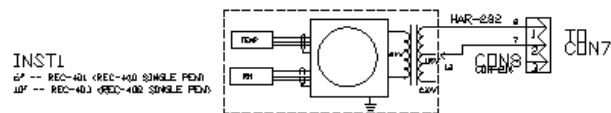


SIDE MOUNT RECORDER


REC-107/103 (RCDR320 & RCDR321)

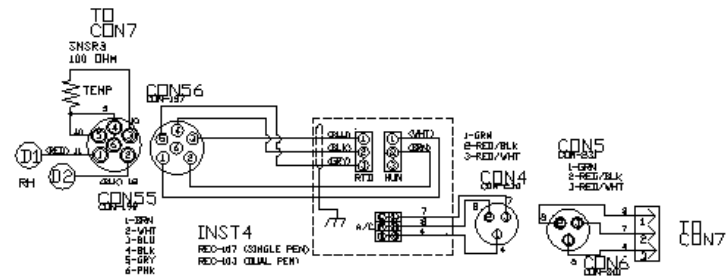

REC-201/202 (RCDR322 & RCDR323)

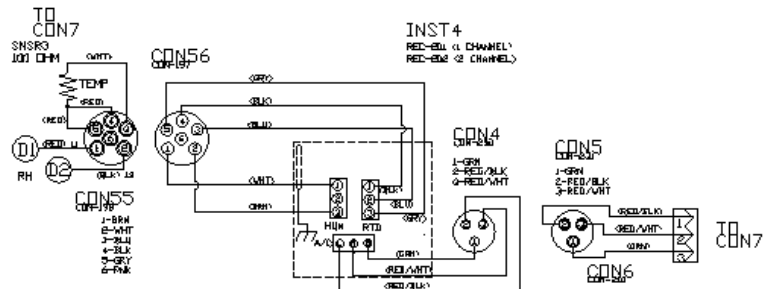

VHP MODULE

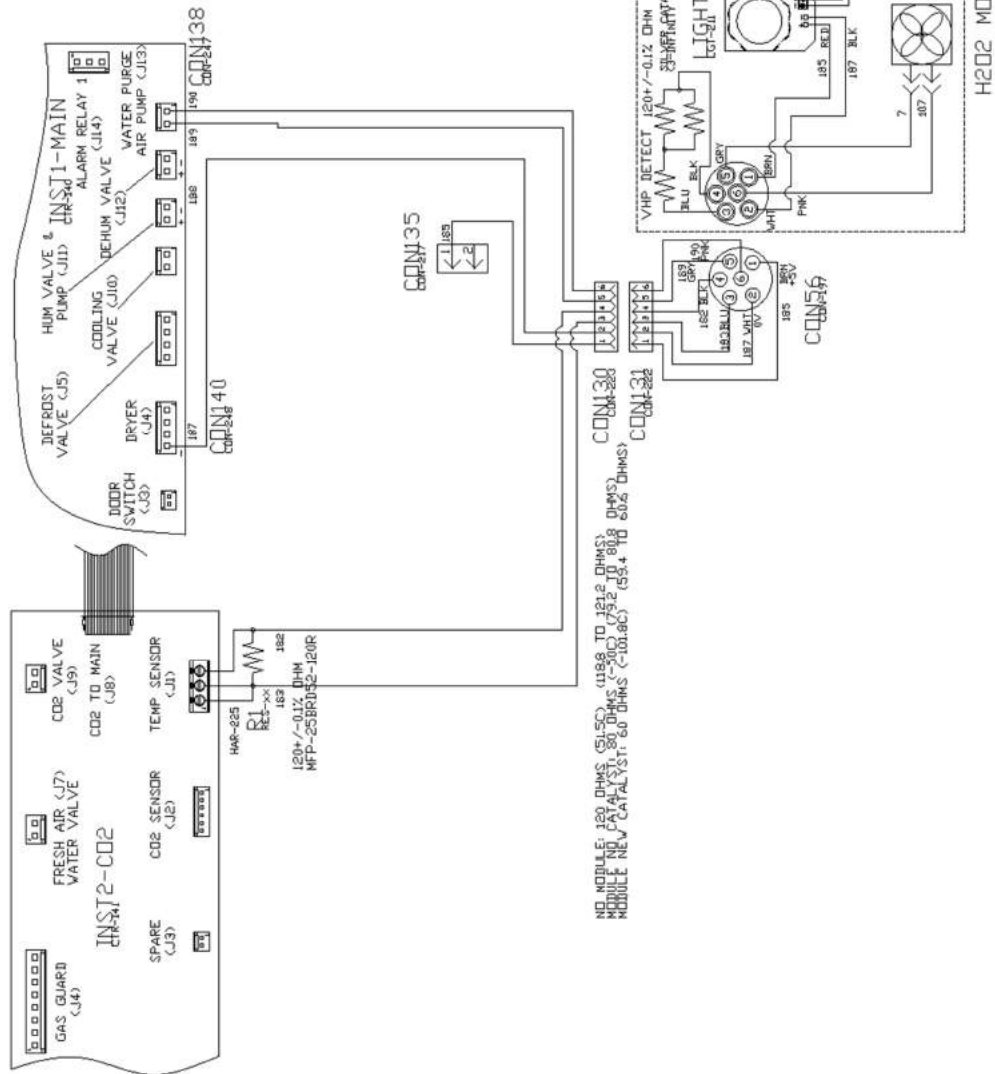

TROUBLESHOOTING

Problem -- Unit will not turn on

- Is the unit connected to a dedicated electrical circuit as defined in the installation section of the manual?
- Is there power at the electric outlet the unit is plugged into?
- Is the unit's power switch turned on?

Problem -- Unit temperature is above / below temperature setpoint

- Has the unit's temperature setpoint been recently lowered / raised and if so has the unit been allowed 12 hours stabilize at the new setpoint?
- Has the door been recently opened for an extended period of time?
- Are the access port stoppers installed in the cabinet?
- Is the condenser filter on the front of the cabinet clean?

Unit humidity level is above / below humidity setpoint (HUMD311-1,-4)

- Is the unit connected to a water source as specified in the installation section of the manual?
- Has the unit been leveled to insure the cabinet drain works correctly?
- The cabinet's drain line uses gravity to remove water. Does the drain line have any rises in it above the cabinet's drain level that could be trapping water?
- Has the unit's humidity setpoint been recently lowered / raised and if so has the unit been allowed time to stabilize at the new setpoint?
- Has the door been recently opened for an extended period of time?
- Are the access port stoppers in the sides of the cabinet installed?
- Is the condenser filter on the front of the cabinet clean?

Lights will not come on?

- Make sure light banks are plugged in to the correct connector
- Check the light timer screen and make sure the touch screen programming is correct

SPARE / REPLACEMENT PARTS



CAUTION: Before servicing the unit, the mains power supply cord must be unplugged to avoid risk of shock. Any area of the unit that requires a tool to access shall only be serviced by trained personnel approved by Caron Products.



R290 REFRIGERANT UNITS

Do not damage the refrigeration circuit. Component parts shall be replaced with like components and servicing shall be done by authorized personnel to reduce the risk of possible ignition.

General

| Part Number | Description |
|-------------|----------------------------------|
| MTR-130 | Blower Motor |
| BLW-112 | Blower Wheel |
| CTR-140 | Main Controller Board |
| POW-115 | 24V DC Power Supply |
| FLTR301 | Condenser Filter Replacement Kit |
| CRD-110 | Power Line Cord (115V 60Hz) |
| CRD-112 | Power Line Cord (230V 60Hz) |
| CRD-108 | Power Line Cord (230V 50Hz) |
| STP-101 | 2" rubber port stopper |



The mains power supply cord must be replaced by the corresponding CRD part number above. The use of an inadequate mains power supply cord could result in equipment failure or personal harm to the user.

Temperature Related

| Part Number | Description |
|-------------|-------------------------------------|
| HTR-153 | Air Heater |
| RMT-117 | 118C Air Heater Thermostat |
| RTD-101 | Temp Sensor -- RTD 100 Ohm Platinum |
| CMP-134 | 115V / 60Hz Condensing Unit |
| CMP-135 | 230V / 50/60Hz Condensing Unit |
| SOL-147 | Refrigeration Cooling Solenoid |

Humidity Related (HUMD311-1,-4)

| Part Number | Description |
|-------------|---------------------------------------|
| FIT-394 | 3/8 Barb Fitting (2X) |
| HUM-110 | RH Sensor |
| PMP-150 | 24VDC RH Pressure Pump |
| NOZ-110 | Precision RH Spray Nozzle |
| SOL-135 | Humidification Solenoid |
| PMP-157 | Air Pump,115V |
| PMP-160 | Air Pump,220V |
| TUB-183 | Drain Tubing, Black, 3/8"ID x 1/2" OD |
| TUB-132 | Water Supply Tubing, Black, 1/4" |

Fuse Related

NOTE: All fuses are slow blow

| ID | Description | 115V | 230V |
|------|-----------------------------|---------------|---------------|
| SW1 | Main circuit breaker switch | CBR-112 (16A) | CBR-115 (10A) |
| FUS1 | Heater fuse | FUS-103 (10A) | FUS-104 (5A) |

* Fuse size varies depending upon whether the chamber has a single internal duplex outlet or two internal duplex outlets

Options Related

| Part Number | Description | Option |
|-------------|-------------------------------------|------------------|
| PEN-103 | Red pen for 6 inch recorder | RCDR316, RCDR317 |
| PEN-104 | Blue pen for 6 inch recorder | RCDR317 |
| PPR-104 | 6 inch recorder paper, 7 day 0-60C | RCDR316 |
| PPR-105 | 6 inch recorder paper, 7 day 0-100C | RCDR317 |
| PPR-106 | 10 inch recorder thermal paper | RCDR318, RCDR319 |
| TUB-174 | 1/2" I.D. silicone tubing | PUMP301 |
| WIR-102 | 20/3 conductor shielded wire | ALRM302 |

Caron Products and Services, Inc.
27640 State Route 7
Marietta, OH 45750 USA

Declares that the product:

Designation: 7340, 7350 Series
Model Numbers: 7340-25, 7340-33-3, 7350-25, 7350-33-3
Classification: Electrical equipment intended for residential, commercial and lighting industrial environments
Rated Voltage: 220-240 ~ (ac)
Rated Frequency: 50Hz

Meets the essential requirements of the following European Union Directive(s) using the relevant section(s) of the normalized standards and related documents shown:

Directives: Low Voltage 2014/30/EU, EMC 2014/30/EU, RoHS 2011/65/EU

Standard: IEC 61010-1:2010

Safety requirements for electrical equipment for measurement, control, and laboratory use. Part 1: General Requirements.

Standard: IEC 61010-2-012:2016

Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 2-012: Particular requirements for climatic and environmental testing and other temperature conditioning equipment

Standard: EN 61326-1:2012

Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements

Signed for and on the behalf of Caron.



Date of issue: April 1, 2021 in Marietta, OH, USA

By: Bob Dotterer
Engineering Manager, CARON

DECLARATION OF CONFORMITY

Caron Products and Services, Inc.
27640 State Route 7
Marietta, OH 45750 USA

Declares that the product:

| | |
|------------------|--|
| Designation: | 7340, 7350 Series |
| Model Numbers: | 7340-25, 7340-33-3, 7350-25, 7350-33-3 |
| Classification: | Electrical equipment intended for residential, commercial and lighting industrial environments |
| Rated Voltage: | 220-240 ~ (ac) |
| Rated Frequency: | 50Hz |

Meets the essential requirements of the following UK legislation using the relevant section(s) of the UK designated standards and related documents shown:

UK legislation: Electrical Equipment (Safety) 2016, EMC 2016, RoHS 2012

Standard: IEC 61010-1:2010

Safety requirements for electrical equipment for measurement, control, and laboratory use. Part 1: General Requirements.

Standard: IEC 61010-2-012:2016

Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 2-012: Particular requirements for climatic and environmental testing and other temperature conditioning equipment

Standard: EN 61326-1:2012

Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements

Signed for and on the behalf of Caron.



Date of issue: December 30, 2021 in Marietta, OH, USA

By: Bob Dotterer
Engineering Manager, CARON